


REV	DESCRIPTION	DRWN	CHKD	APPR
001	MSM Series 100 & Series 200 Limits	TC	IM	TC

NOTES:

Positions of Standard Petri Dish Limits for Default "Singer" Layout MSM Series 100 & 200
See Section 5.6 Limits in User Handbook

Note: when Changing Limits, work Clockwise starting at NE 1 Drive stage using joystick until Display shows correct values, (the display will update on release of the joystick, so trial and error until desired position is shown then press the Equals Key to store that value, then continue procedure until all eight Limit points are set. When finished, press Arrow key to exit the Limits Loop.


MATERIAL: XXX		FINISH: XXX	
PART NAME: MSM 200 Plate Limits			
PART NUMBER: XXX		SHEET 1 - 1	©
DRWN	DATE	CHKD	DATE
TC	02/10/14	IM	02/10/14
APPR	DATE	APPR	DATE
TC	02/10/14	TC	02/10/14



**SINGER
INSTRUMENTS**
A RESPONSIBILITY TO SCIENCE!
ROADWATER WATCHET SOMERSET TA23 0RE UNITED KINGDOM
TEL 00 44 1984640226

GLOBAL TOLERANCES UNLESS OTHERWISE STATED TO BE TO SINGER GENERAL TOLERANCE SHEET

ALL DIMENSIONS IN MILLIMETERS

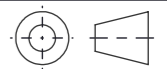

DO NOT SCALE
IF IN DOUBT ASK