

ROTOR – DELETING MANUAL ROTINES


INTRODUCTION

The ROTOR Manual operation allows you to create routines (programs) which makes the ROTOR work the way you want it to. It is a very easy system to use and it will have been demonstrated during the commissioning of your ROTOR.


This document explains how to delete the routines you have created.

PROCEDURE

- Select the “i” (information) on the right side of the screen.

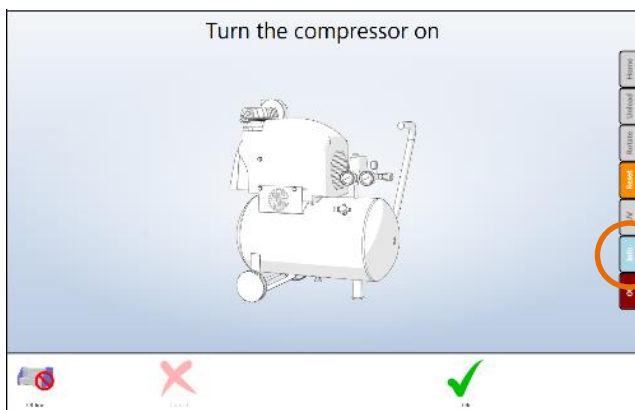


Laptop MCI

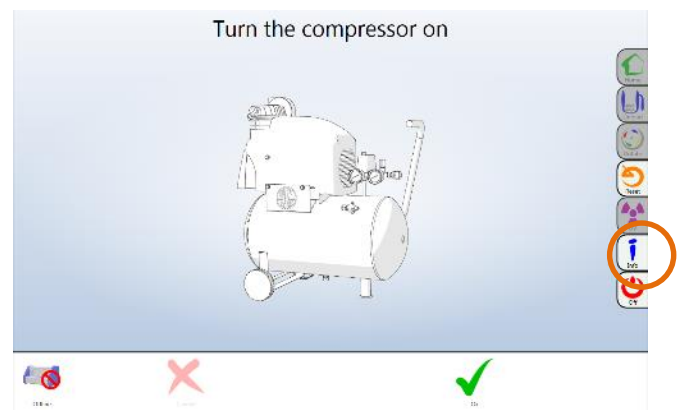

Touch screen MCI


- Select “Manage programs in your library”. This facility is also available by selecting the ‘Advanced Programming’ option in the programming options menu.

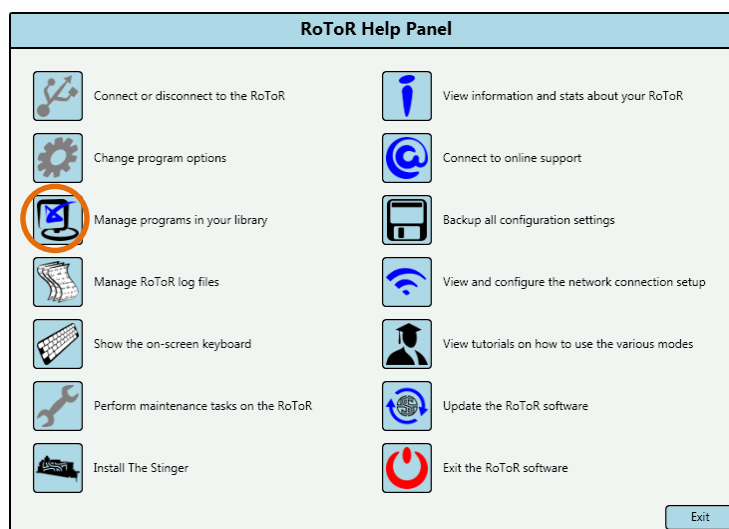

Manage programs

ROTOR – DELETING MANUAL ROTINES

- The screen shown below will be displayed. The tab “Delete” will be highlighted and a list of all your routines will displayed. Select the box next to the routine you wish to remove or the one at the bottom of the screen “select all programs”. When the file you wish to delete is selected there will be an ‘X’ in the box.

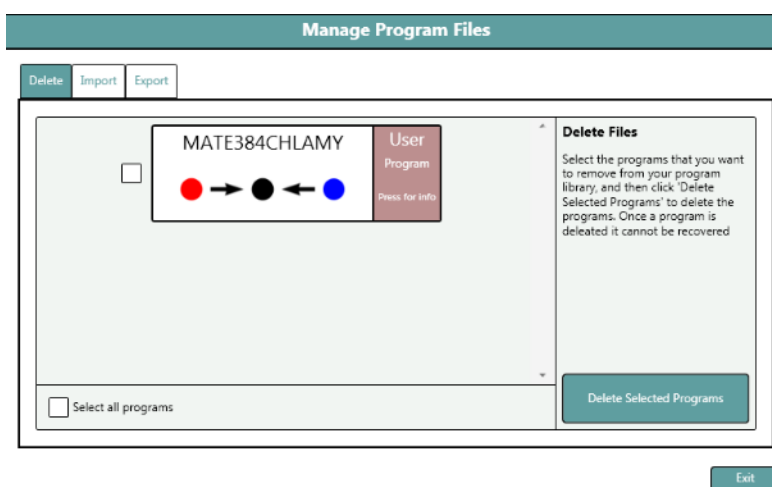

- Once you are happy with your selection simply select the button “Delete Selected Programs”.
- Select “Exit” and you will get a message confirming the routines have been deleted. This will be followed by a message asking if you wish to re-buffer the SRR routines to which you answer “OK”.
- You can then exit out of the Help Panel.
- The routines have been deleted.

CONCLUSION

We hope that this explanation will assist you in performing the above procedure. If you need further assistance do not hesitate to contact the Technical Support Team at Singer Instruments.

CONTACT INFORMATION

Name(s): *Ian Morgan*

Phone : *+44 (0) 1984 640226*

Email: *technicalsupport@singerinstruments.com*