

PACKING THE SINGER MSM SYSTEM 300

INTRODUCTION

The following is a brief instruction for packing up your MSM System 300 for transport. If you still have the original transit brackets we would advise you to fit them.

PROCEDURE - Switch off the instrument and isolate from the mains.

- **All cables have to be removed from the MSM**
 - a) Umbilical (25 way) cable and connectors.
 - b) Joystick control cable and connectors.
 - c) Micro-Zapper, Mini-Din Lead.
- **Casio PDA**
 - a) Disconnect Casio 'D' Type connector from the MSM.
 - b) Remove the connector from the PDA.
 - c) Remove the power supply lead from the PDA.
- **Trinocular Head**
 - a) Take eyepieces out of their sockets.
 - b) Unscrew the locking clamp until slight resistance is felt.
 - c) Remove trinocular head by pushing it to the right against the sprung loaded catch.
- **Objectives**
 - a) Unscrew x20 objective and place into plastic receptacle with screw cap.
 - b) Unscrew x4 objective and place into plastic receptacle with screw cap.

(If you no longer have the plastic receptacles use small plastic bags)
- **Micromanipulator**
 - a) Release the lever by turning it anti-clockwise while holding the weight of the micromanipulator. Slide it out of the dovetail and place it with the other items removed.
- **Joystick Control**
 - a) Release this in the same way as the micromanipulator and place it with the other items.

CONCLUSION

We hope that this explanation will assist you in performing the above procedure. If you need further assistance do not hesitate to contact the Technical Support Team at Singer Instruments.

CONTACT INFORMATION

Name(s): *Ian Morgan*

Phone : *+44 (0) 1984 640226*

Email: *technicalsupport@singerinstruments.com*