




SINGER
INSTRUMENTS
A RESPONSIBILITY TO SCIENCE!

PHEN~~Ø~~BOOTH+™

USER GUIDE v1.4

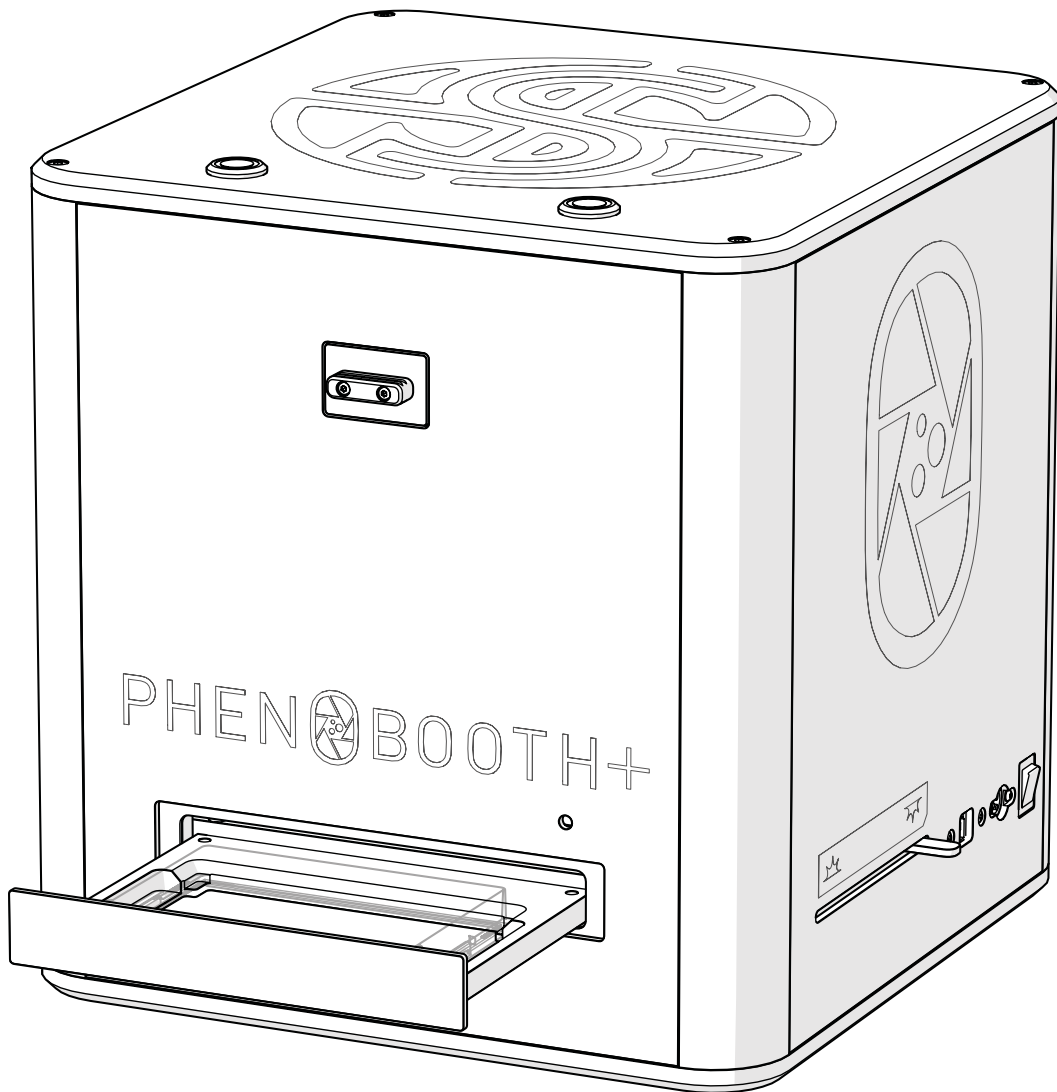



READ BEFORE YOU START!

- **Remove plate lids before imaging.** If lids are in place when the input tray is closed, the machine will crash and may be damaged.
- **Keep your hands clear while the input tray is moving.** Catching your hands in the input tray may cause injury.
- **Ensure power supply is the correct way round.** Plugging the power supply into the PhenoBooth+ upsidedown will cause damage to the machine. See p.10 for details.
- **Avoid plugging the USB 3.0 cable in slowly.** A common error with all USB 3.0 devices is that they are recognised as USB 2.0. This happens when the USB plug is inserted too slowly.
- **Remove plate and petri adaptor before moving PhenoBooth+.** If you need to move the PhenoBooth+ for any reason, make sure any plates are removed from the machine as well as the petri dish adaptor. These items can fall from the input tray and become lodged inside the PhenoBooth+.
- **The PhenoBooth+ is not a microscope!** It is a macroscopic CCD imaging workstation. As such it will not detect sub-cellular or low levels of expression. The in-house strains used to successfully test PhenoBooth's fluorescence detection have a fluorophore molarity of 1.84×10^{-30} nM per cell. As such we cannot guarantee fluorescence detection below this value.
- **Set agar height for different plate types.** The PhenoBooth+ is capable of using a variety of different plates which can have different base thicknesses. To account for this, we ask users to input the distance from the top of the agar (trying to ignore the meniscus) to the bottom of the plate. See p.53 for details.

2.	Before you start!	32.	RCP PROJECT
3.	Introduction	33.	New project
5.	OUT OF THE BOX	34.	Image acquisition
6.	Anatomy & features	36.	Pre-processing & detection
9.	Unboxing & installation	37.	Filtering
11.	LET'S START!	38.	Stinger Export
12.	PhenoBooth+ hardware operation	39.	ROTOR HDA / The Stinger
	PhenoBooth+ software set-up	43.	BEST PRACTICE
14.	COLONY COUNTING PROJECT	44.	Software icons
15.	New project	46.	Feedback light
16.	Image acquisition	47.	Batch processing
17.	Pre-processing	48.	History plug-in
19.	Detection	49.	Add new plate run
20.	Filtering	50.	Import/export Project
21.	Export	52.	Using filters
22.	COLONY SCREENING PROJECT	53.	Setting Agar Height
23.	New project	54.	Notes
24.	Image acquisition		
26.	Pre-processing		
27.	Detection		
28.	Filtering		
29.	Analysis		
30.	Visualisation		
31.	Export		

PHENOBOOTH+™

INTRODUCTION

The PhenoBooth+ is a fast, semi-automated plate reader for imaging and analysis of cell colonies. The PhenoBooth+ can image standard SBS plates, Petri dishes and Singer PlusPlates.

USER GUIDE

Follow these instructions to get the most out of the PhenoBooth+. This guide outlines getting started and general operation, as well as software tutorials for each package. The information in this guide relates to software version 2.22.0607.1

DISCLAIMER


At Singer Instruments, we are constantly seeking to improve our products and adapt them to the requirements of modern research techniques and testing methods. This involves modification to the mechanical structure and optical design of our instruments. Therefore, all descriptions and illustrations in this user guide, including all specifications are subject to change without notice. To download the latest version of this user guide please visit: singerinstruments.zendesk.com

OUT OF THE BOX

Find out what comes with the PhenoBooth+. We'll take you through the steps involved in unboxing and assembling ready for the exciting world of colony imaging!

ANATOMY & FEATURES



PHENOBOOTH+

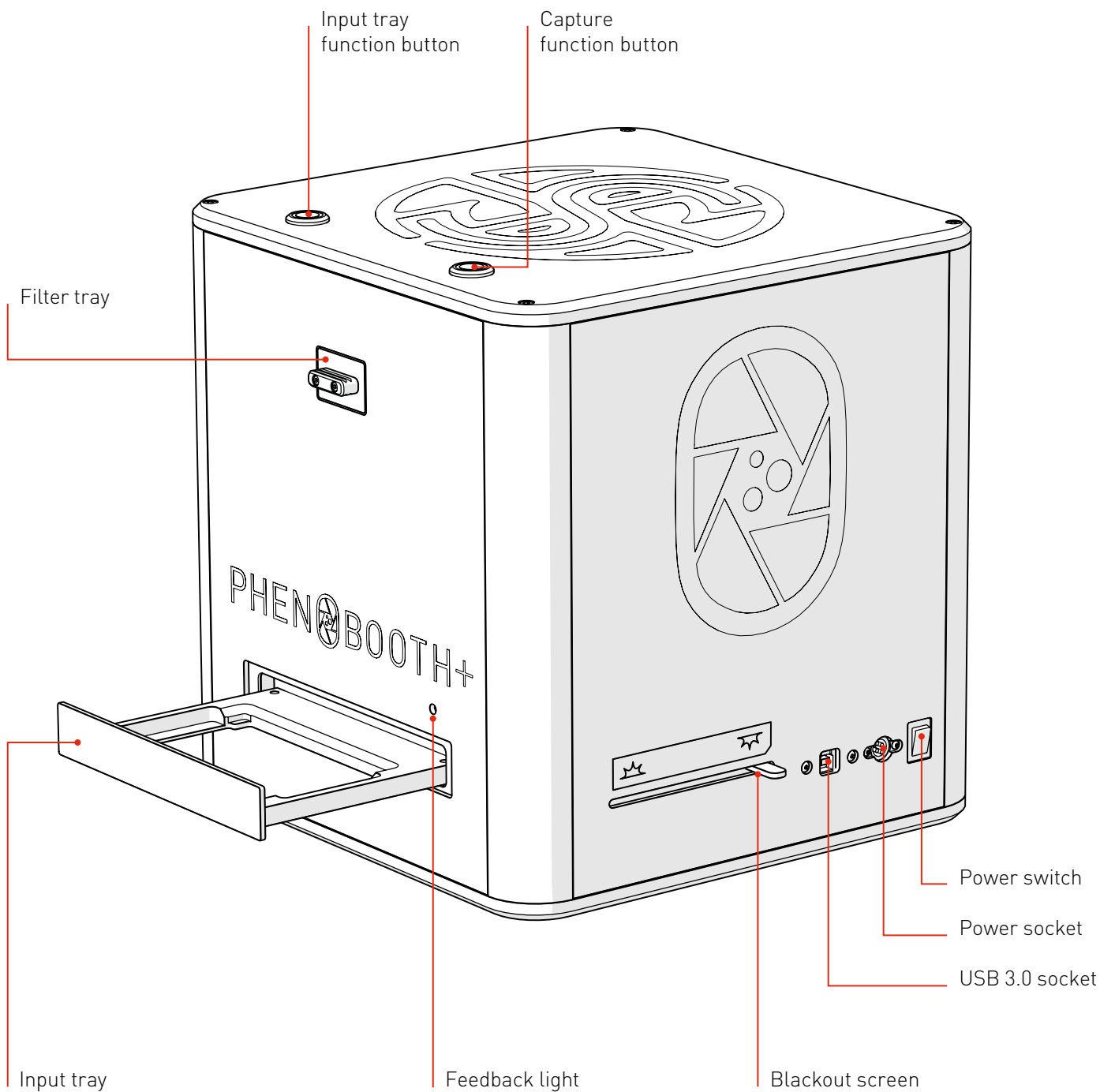

ANATOMY & FEATURES


COMPUTER AND MONITOR

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See Dell instructions in product boxes.

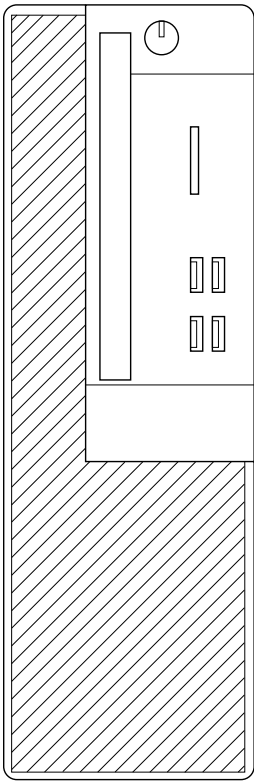

ANATOMY & FEATURES




SOFTWARE

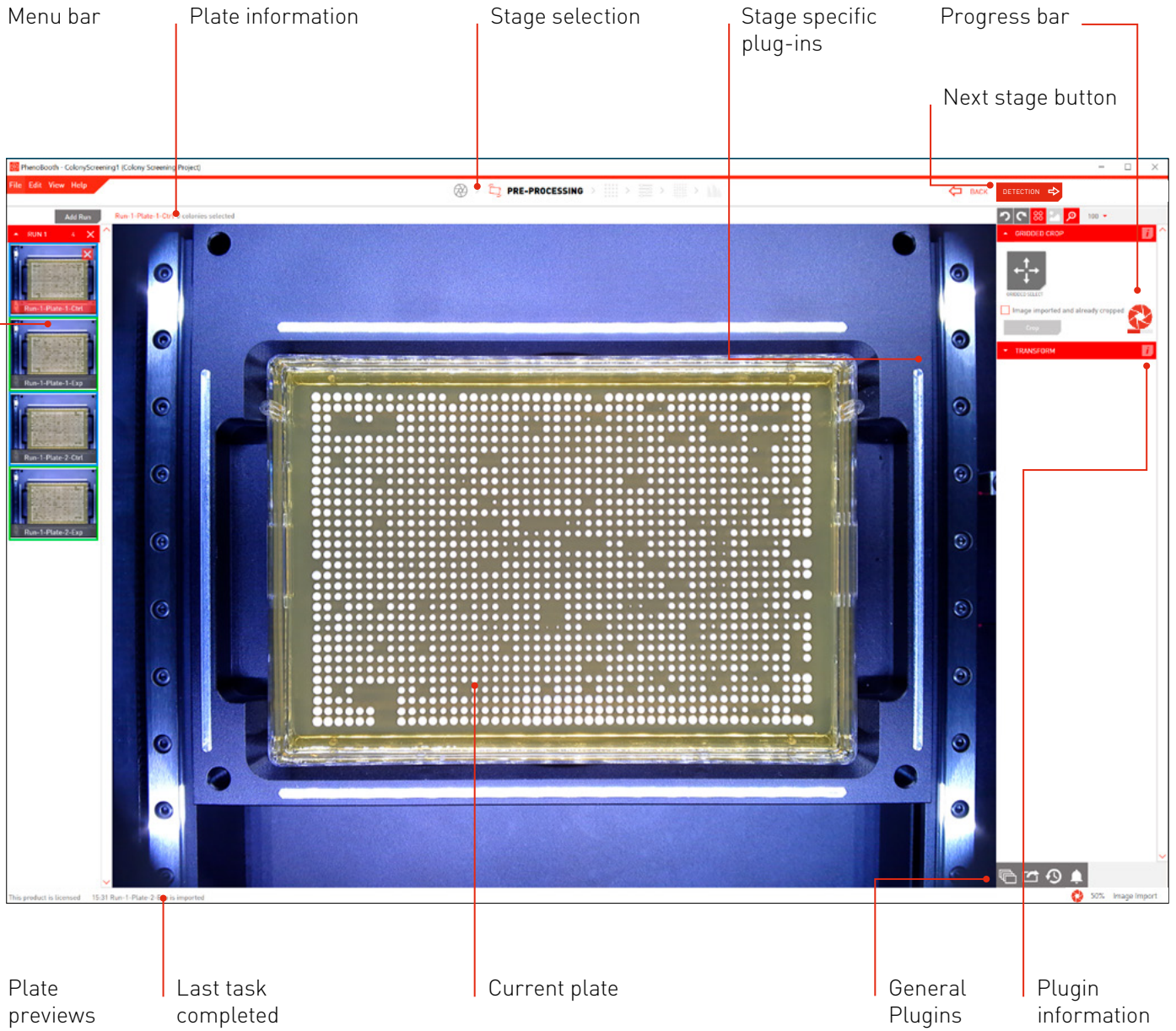

UNBOXING & INSTALLATION

PHENOBOOTH+ BOX

1. PhenoBooth+
2. Power supply
3. 90mm petri adapter (optional extra)

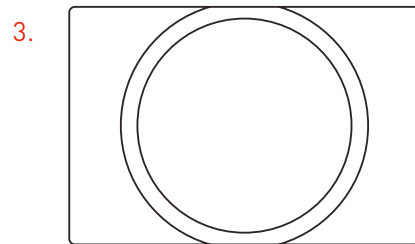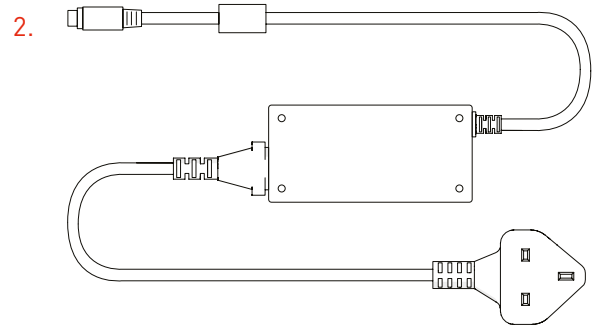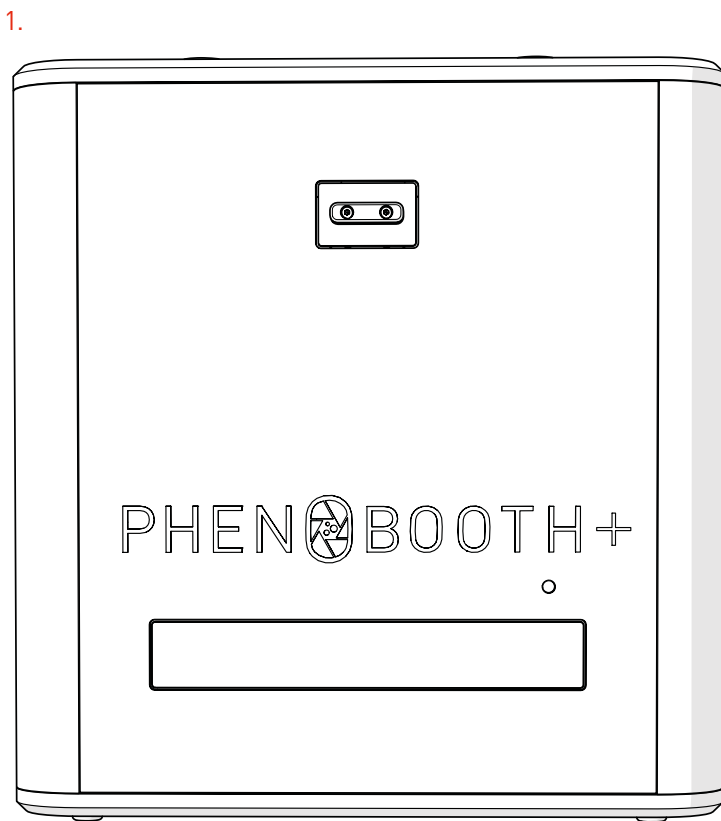


ADDITIONAL ITEMS

4. Monitor box (includes Dell monitor, stand and connections)
5. Computer box (include Dell computer, power cable, mouse and keyboard)


UNBOXING & INSTALLATION

1

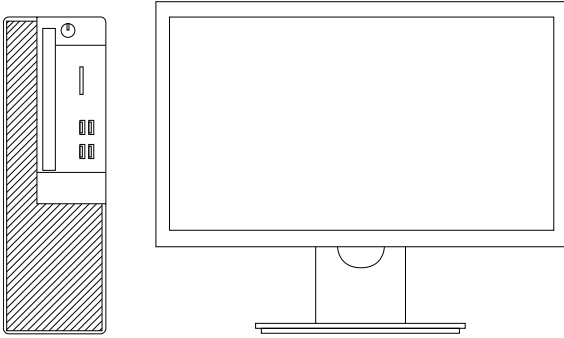


- Set up the computer and monitor as details in their instructions.

3

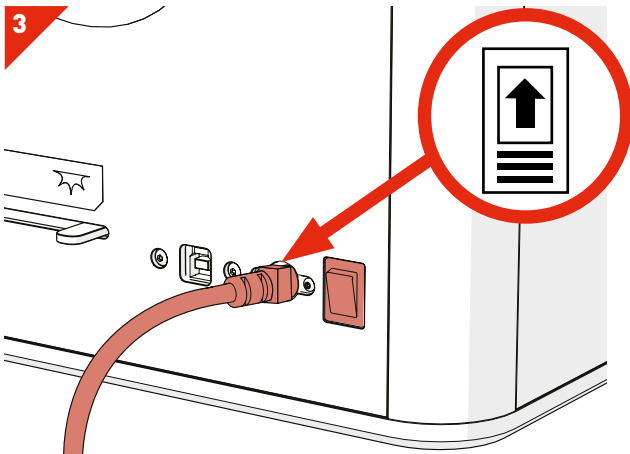


- Insert PhenoBooth+ power supply and turn on the power switch.
- **WARNING:** Make sure the arrow diagram is on top of the cable. Inserting the power supply upside down will cause damage to the PhenoBooth+.

5

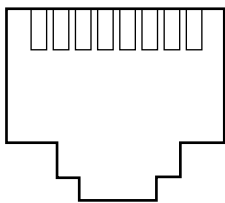


- For future software updates you will need to connect the computer to the internet via the LAN socket. See computer instructions.

2

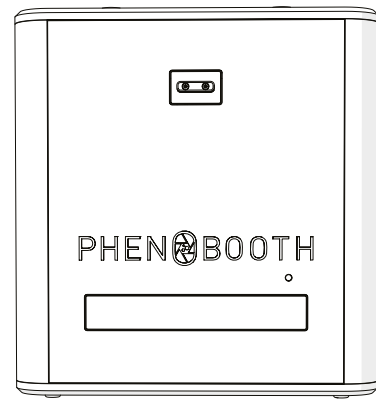

- Carefully remove the PhenoBooth+ from the box and stand on your workspace.
- **HEAVY:** May require 2 people.

4

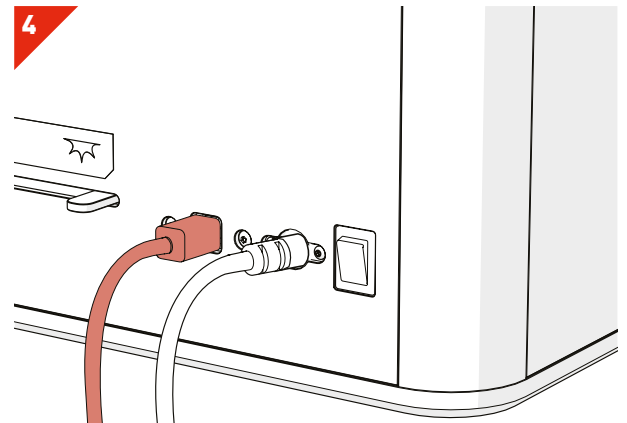


- Insert the USB cable into the PhenoBooth+ USB 3.0 socket and to the MCI USB 3.0 socket.

LET'S START!


I'm ready to start!

Here is a quick run-through of how to use the PhenoBooth+ and get started with the software — it's all very simple!


PHENOBOOTH+ HARDWARE OPERATION

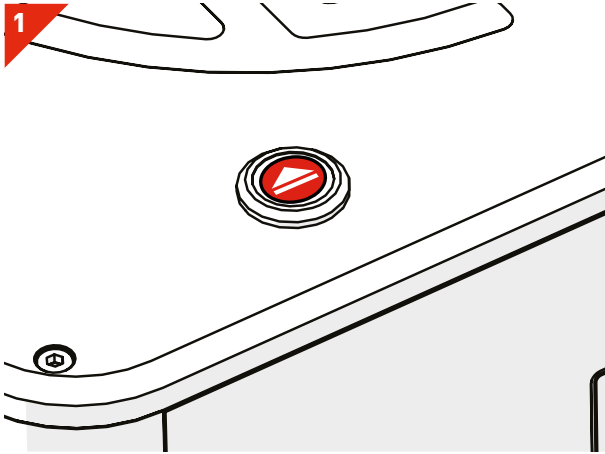


- Press the input tray function button to open the input tray.

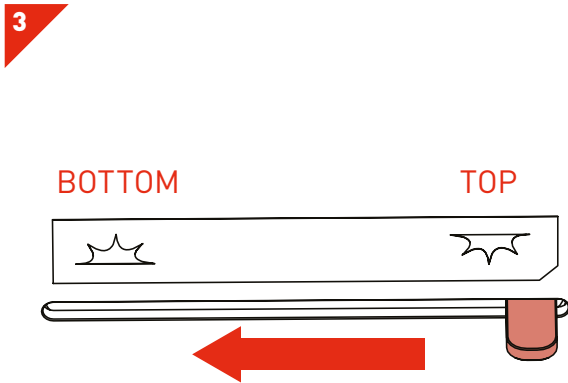


- Choose between top or bottom lighting by sliding the blackout screen (left for bottom and right for top).

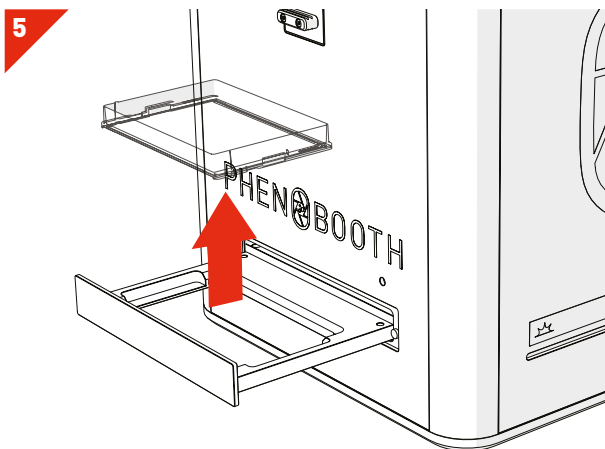


- Press the input tray function button to open the input tray, remove your plate and replace the lid.
- Repeat these steps for the rest of your plates.

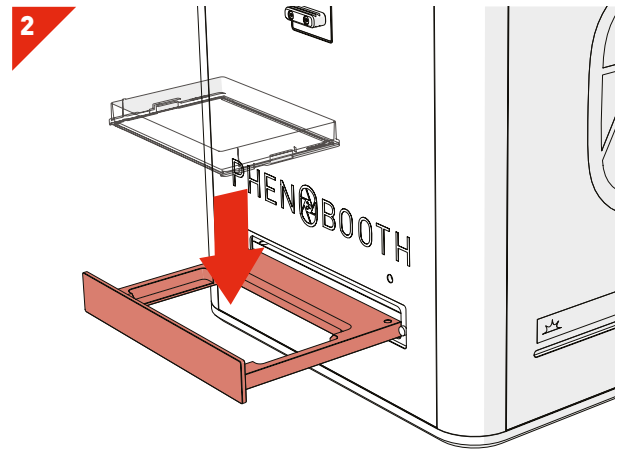

- Remove the lid from your plate and place the plate into the input tray.
- **WARNING:** Ensure plate is nestled correctly and lid is removed to avoid a hardware crash.

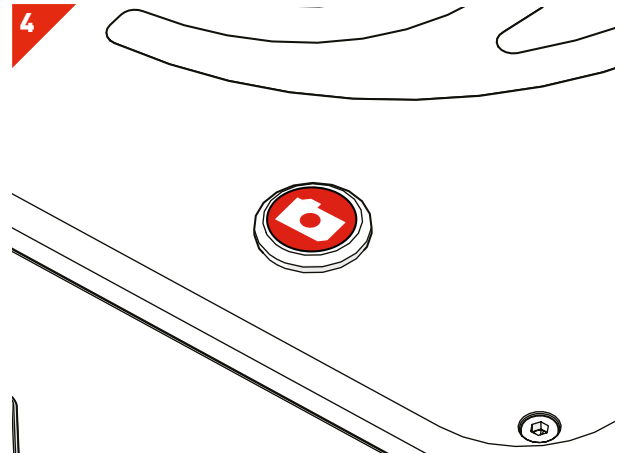


- Press the capture function button to image the plate.
- Follow the software instructions.

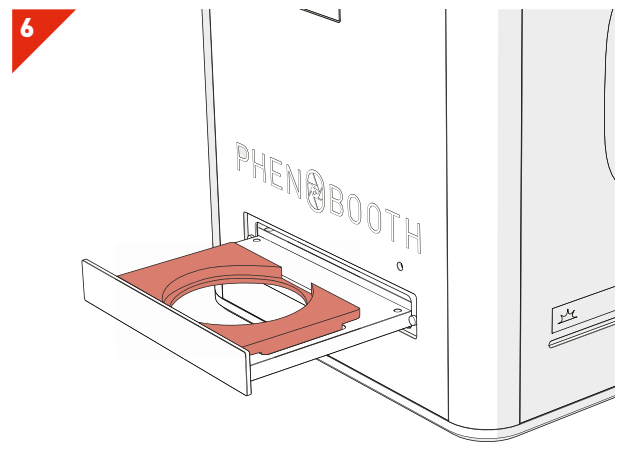

- If using petri dishes, you will need to use a petri dish adaptor. Place the adaptor into the input tray.
- Remove the petri dish lid and insert into the adaptor.


PHENOBOOTH+ SOFTWARE SET-UP

1

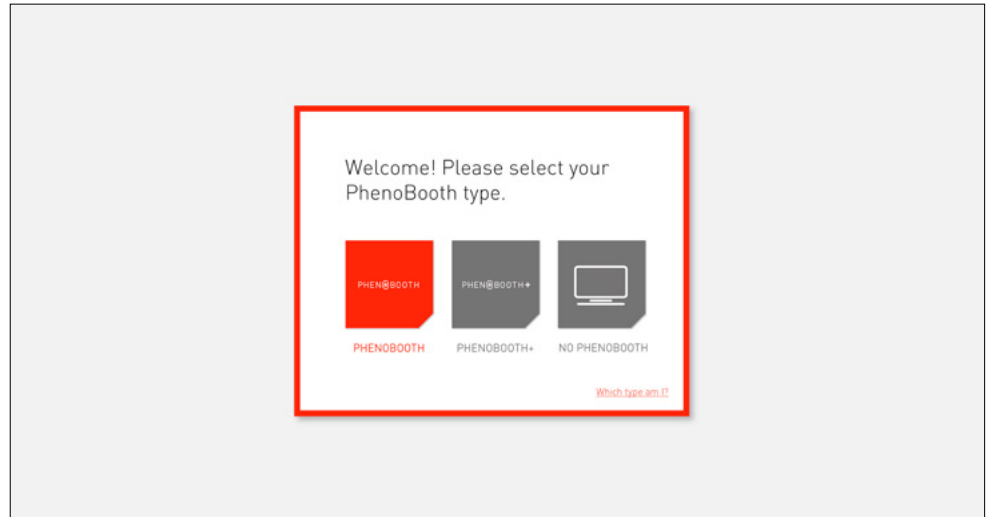


- PhenoBooth+ replaced the original PhenoBooth in October 2019. In order for your software to work properly, you'll need to select your PhenoBooth type:
- **PhenoBooth:** For machines delivered after October 2019
- **PhenoBooth+:** For machines delivered after October 2019. If you're unsure, the label on the front of the machine will tell you which type you have.
- **No PhenoBooth:** For software installed on a stand alone computer.

2

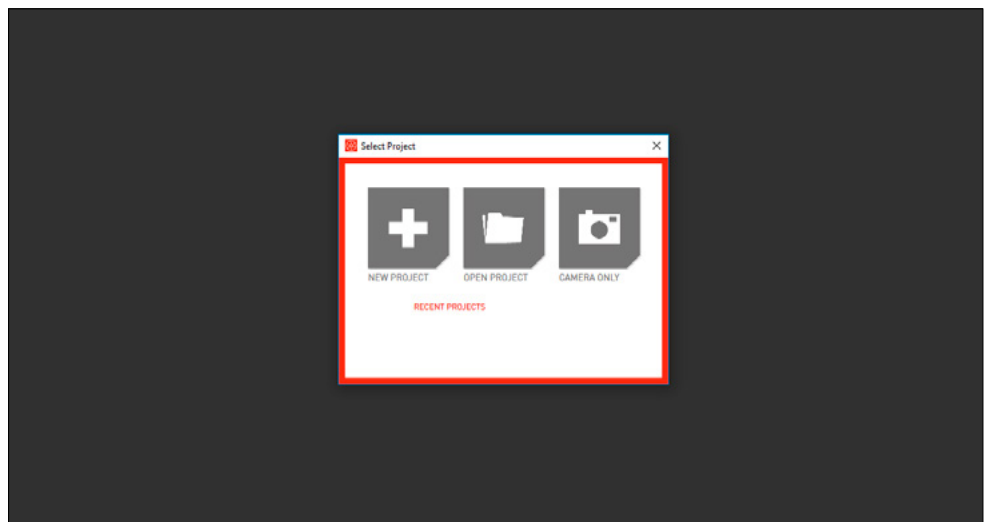


- Now you're all ready to start. Simply click to select a new project or open an existing one.
- If you just want to take a photo of a plate, with no analysis, select Camera Only.

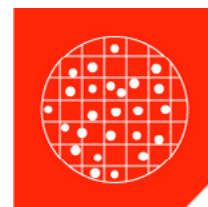

COLONY COUNTING PROJECT



Colony Counting is the basic package included with your PhenoBooth+. There are 5 stages in the Colony Counting package: Image Acquisition, Pre-processing, Detection, Filtering and Summary.

Start a new project then select Colony Counting.

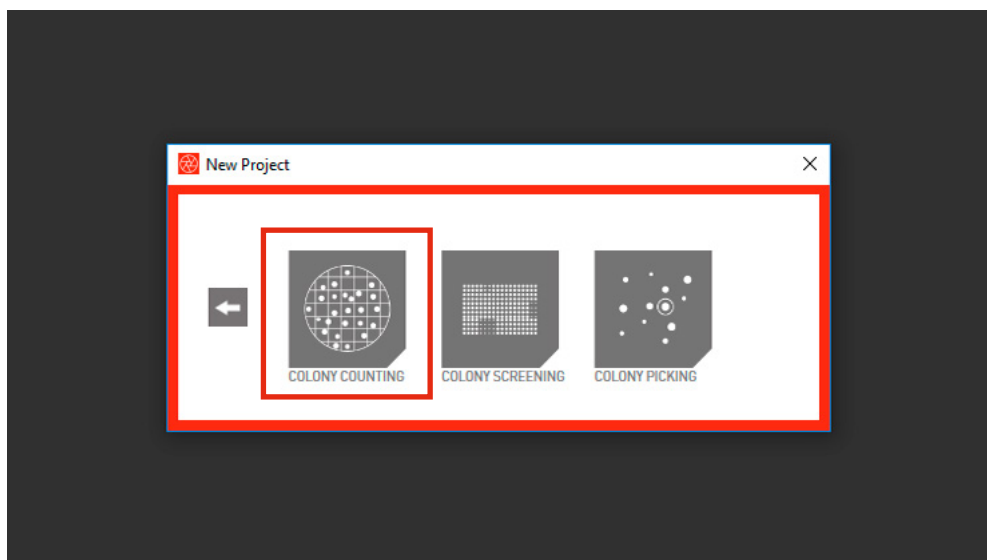



COLONY COUNTING - NEW PROJECT

1

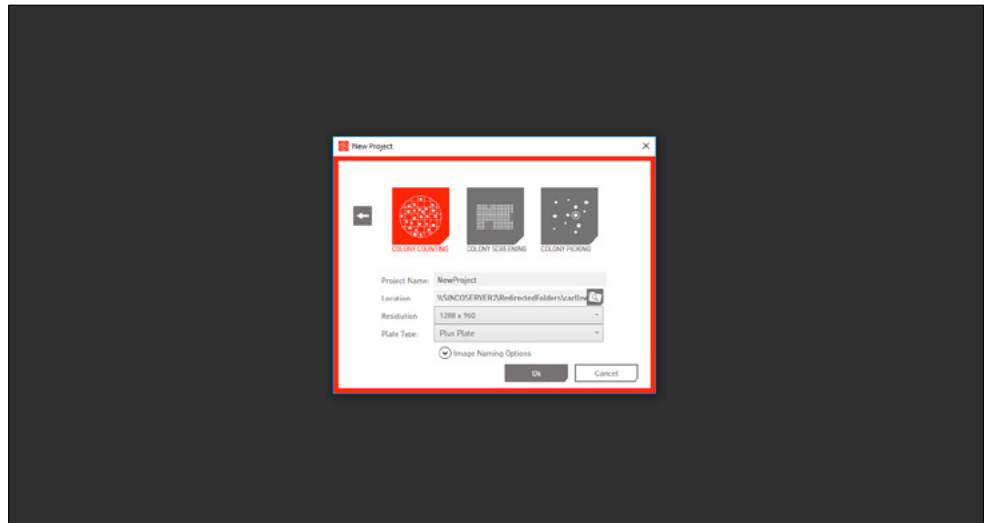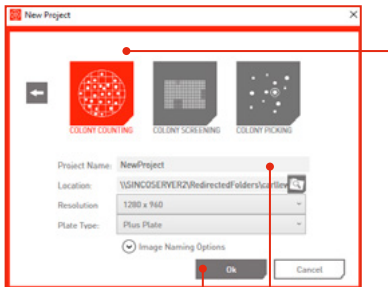

- Open the PhenoBooth+ software and select New Project.
- Select the Colony Counting Icon. Icons will vary depending on which software packages you have purchased. Your chosen project type is highlighted in red.
- Name your project, select a save location, choose an image resolution and select your plate type.
- There are also a few optional naming settings you can use to customise your plate images.
- Click OK to launch project.



2

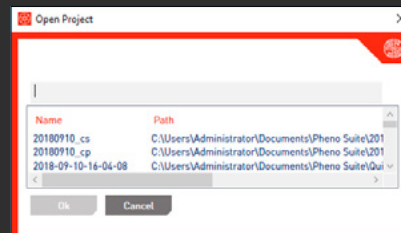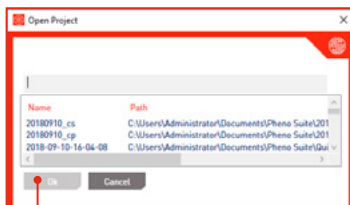

- If you want to open a previous project select Open Project and pick from the list of existing projects.
- Click OK to launch project.


COLONY COUNTING - IMAGE ACQUISITION

1

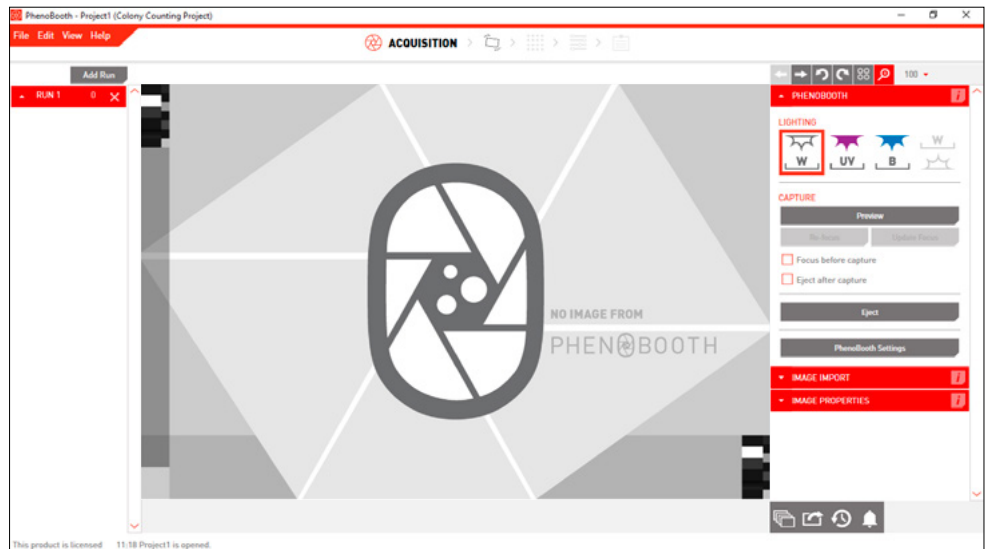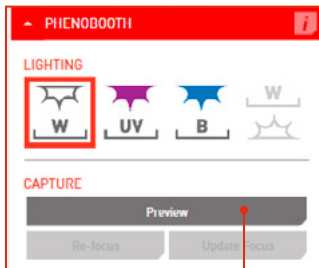

- There are two options for acquiring an image: PhenoBooth+ or import an existing image.
- Select the desired method from the plug-in panel on the right of the screen.
- If using the PhenoBooth+ camera, you can preview the image by pressing Preview.


2

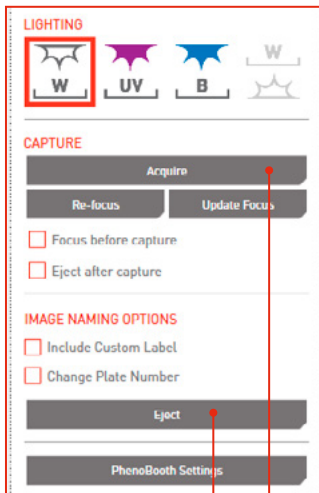

- If using the PhenoBooth+, click on that tab and select your desired light settings. Plate names are automatically generated.
- Insert a plate into PhenoBooth+ and press Acquire.
- Eject will allow you to remove the plate.
- Repeated until all of your plates have been captured.
- ➡ moves onto the pre-processing stage.


COLONY COUNTING - IMAGE ACQUISITION

3

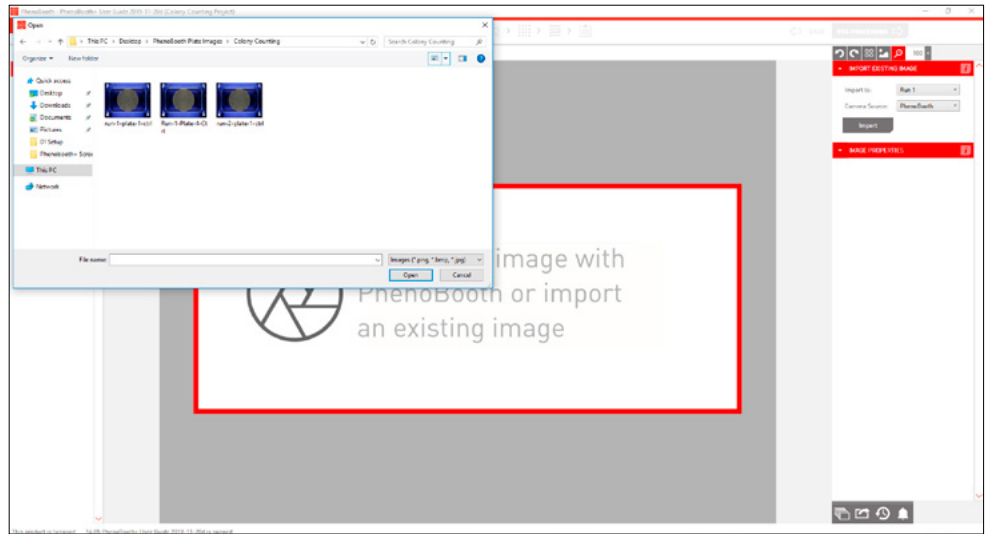

- If using existing images, click the Image Import tab. Here you can select which run the plate images will be associated. If this is your first run then Run 1 will be the only option. See p.49 to add new runs.
- Click Import to select images from your computer.



4

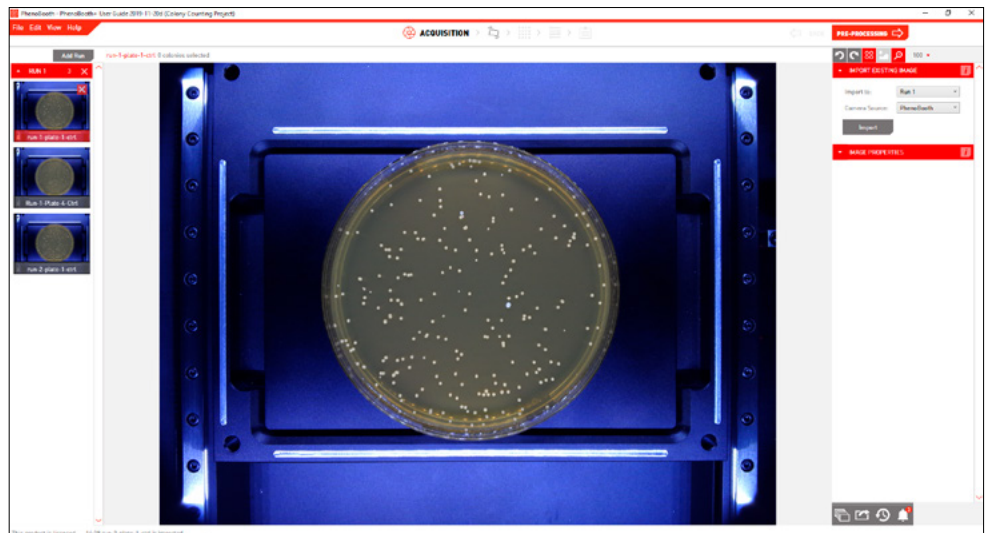

- Your captured or imported images now appear on the left of the screen. The currently selected plate is highlighted with red.
- Once you've finished capturing or importing images, use the next stage button to move onto the pre-processing stage.



COLONY COUNTING - PRE-PROCESSING

5

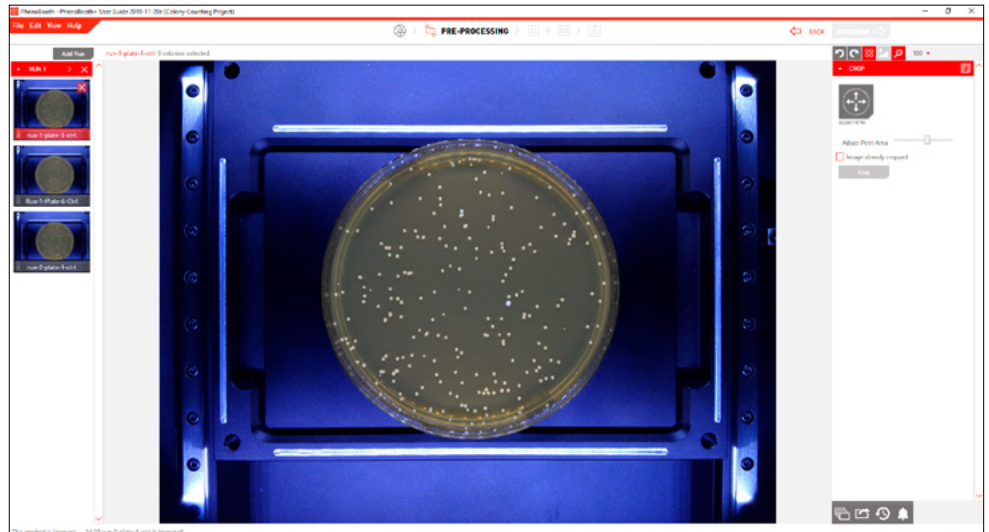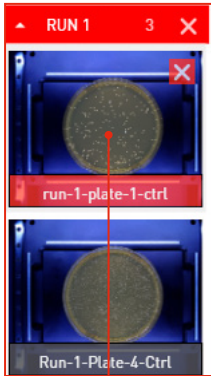

- Select which image to process first. If using the batch process option, this plate's settings will be applied to the rest of the plates in the run. Learn more about batch processing on p.47.



6

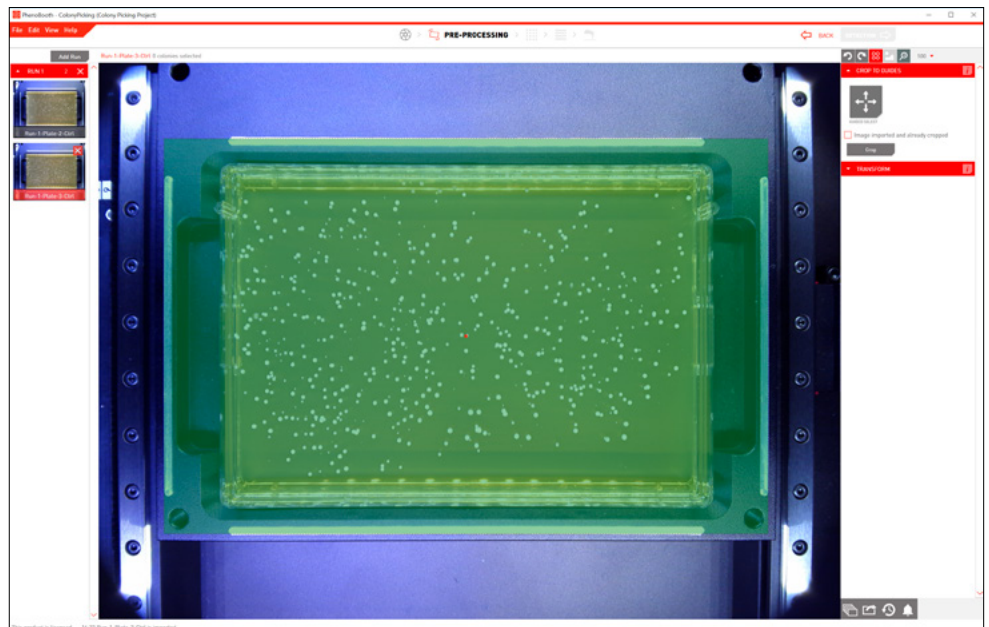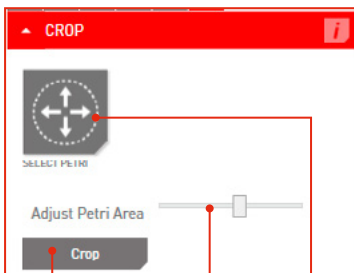

- Click on Select Petri to automatically detect the plate area. You can adjust the plate area using the Petri Slider.
- If using a PlusPlate, press the Select button and drag over the colony area of the plate.
- Click Crop to remove the edge of the plate and beyond.
- You will automatically move onto the filtering stage.



COLONY COUNTING - DETECTION


7

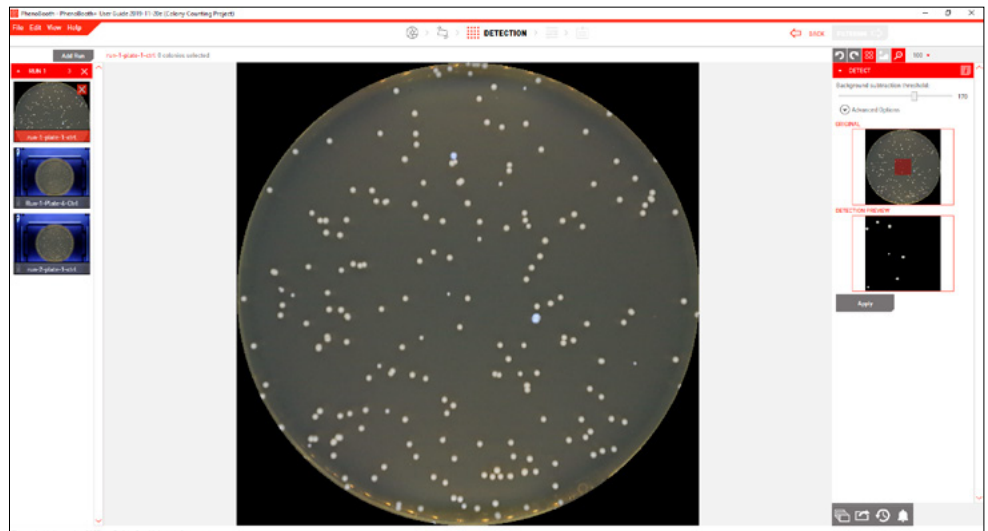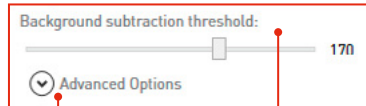


- In order to generate colony information, the background needs to be removed.
- Use the slider to adjust the filter intensity if necessary.
- If you have colonies that touch you can use the colony separation tool found in the advanced options.
- Check that all the smaller colonies are visible in the preview window then click Apply.
- Once applied, you will automatically move onto the filtering stage.

8

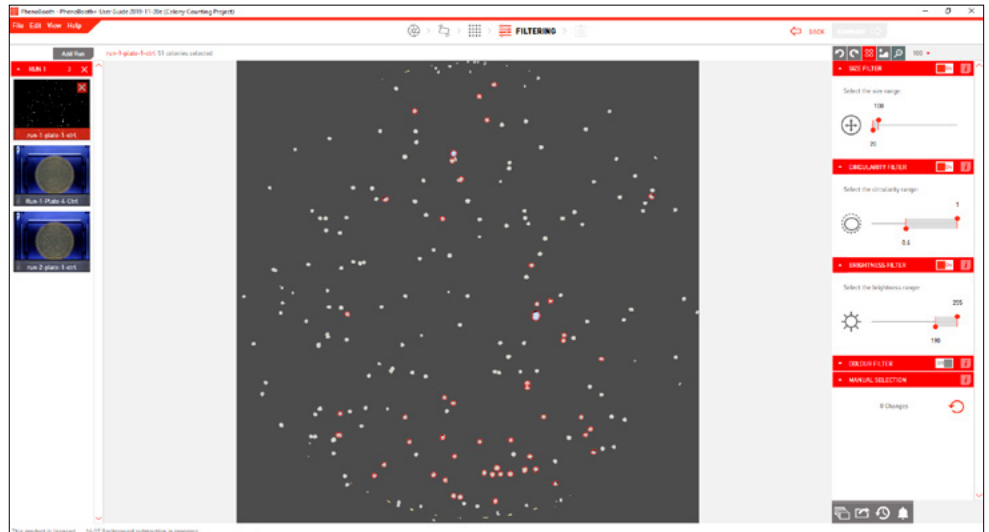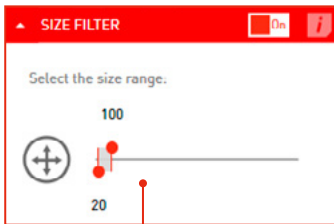

- Detected colonies are highlighted with a red circle.  turns highlights on or off.
- Use the filter options to set minimum and maximum values for size, circularity, brightness and colour for colony detection. You can use the size filter to discard items that are too large or small to be colonies.
- Colonies can also be manually selected by simply clicking on a colony to select or deselect it.




9

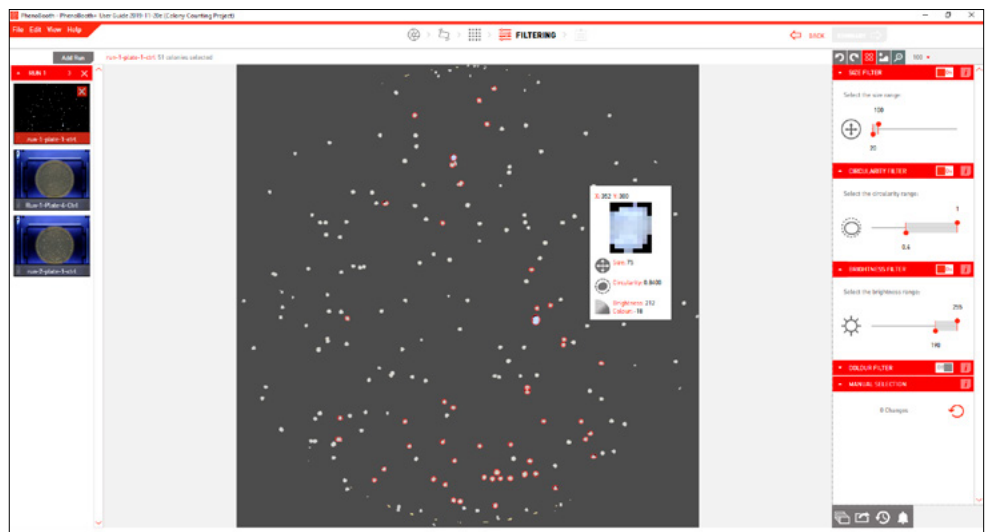

- Hover over a colony to see it's size, circularity, brightness and colour properties in the Colony Info card.
- You can batch process the rest of your plates with the same options using the instructions on p.47.
- After processing all of your plates you can use the next stage button to move onto the summary stage.
- **NOTE:** All plates must be processed before you can advance to the summary.

COLONY COUNTING - SUMMARY

10

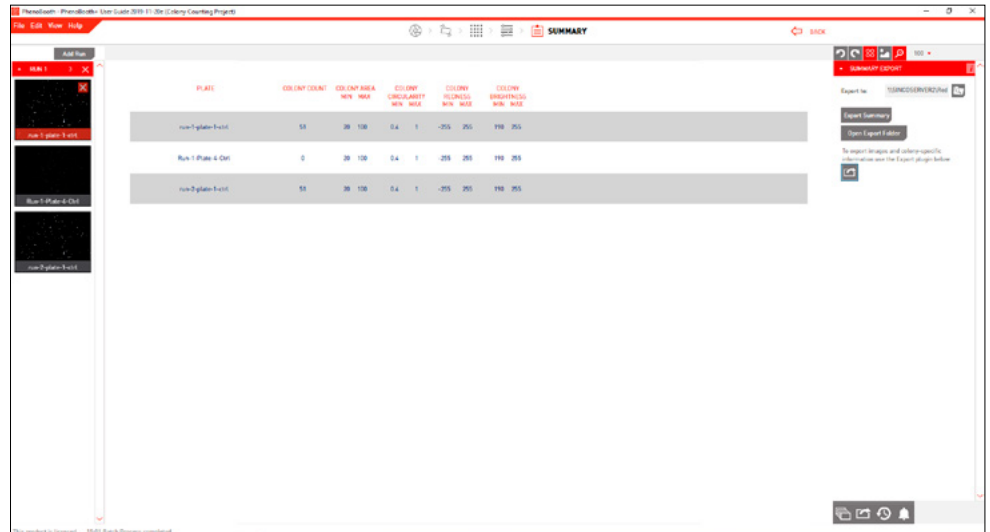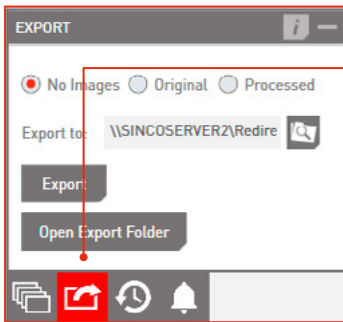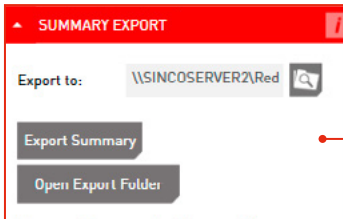

- The Summary Export plugin will export the data on the summary screen.
- For a full export of your workflow, click the Export icon on the bottom right. You can choose to export no images, processed images or the original images.
- The export includes: plate image (if selected), summary file and colony data table.
- The Summary File contains colony counts for all plates as well as colony imaging and filtering options information.
- The Colony Data Table contains position and properties of all detected colonies on all plates.




COLONY SCREENING PROJECT

The Colony Screening package includes all the plug-ins you need for colony screening, such as size and circularity filters, plate normalisation, gene mapping, control / experiment comparison and heat map visualisation.



There are 6 stages in the Colony Screening package: Image Acquisition, Pre-processing, Detection, Filtering, Analysis and Visualisation.

Start a new project then select Colony Screening.

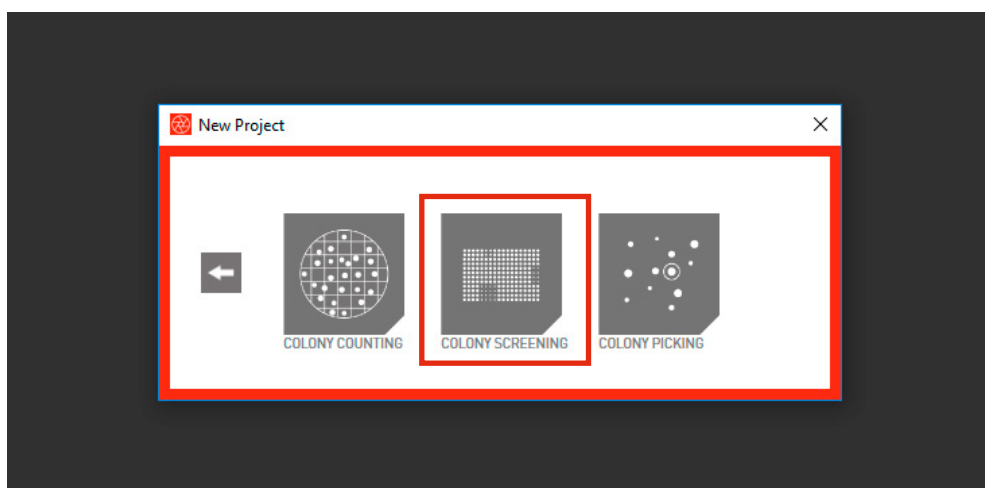



COLONY COUNTING - NEW PROJECT

1

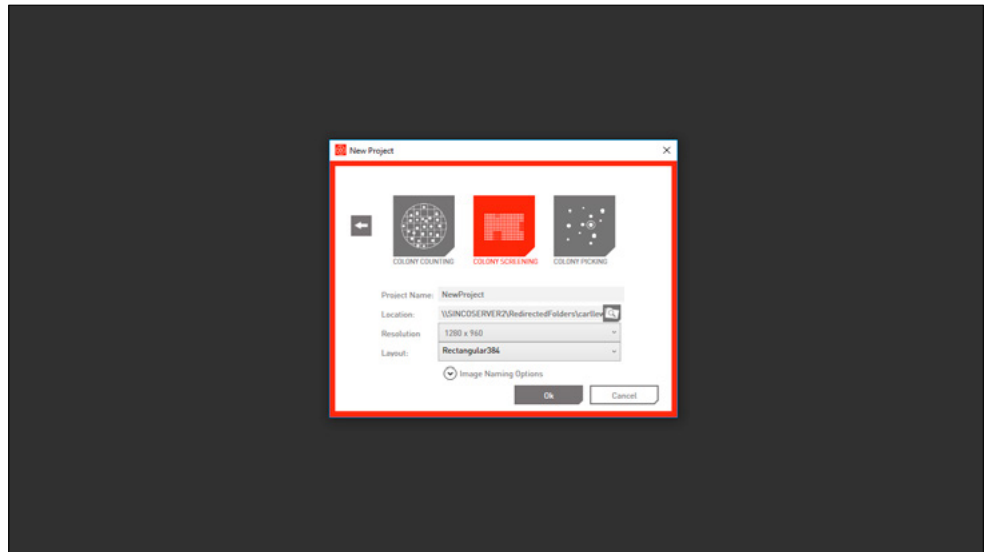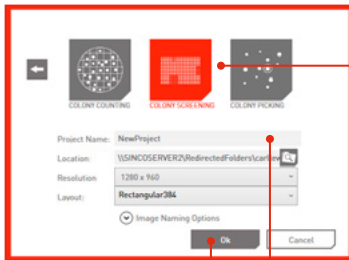

- Open the PhenoBooth+ software and select New Project.
- Select the Colony Screening Icon. Icons will vary depending on which software packages you have purchased. Your chosen project type is highlighted in red.
- Name your project, select a save location, choose an image resolution and your layout.
- There are also a few optional naming settings you can use to customise your plate images.
- Click OK to launch project.



2

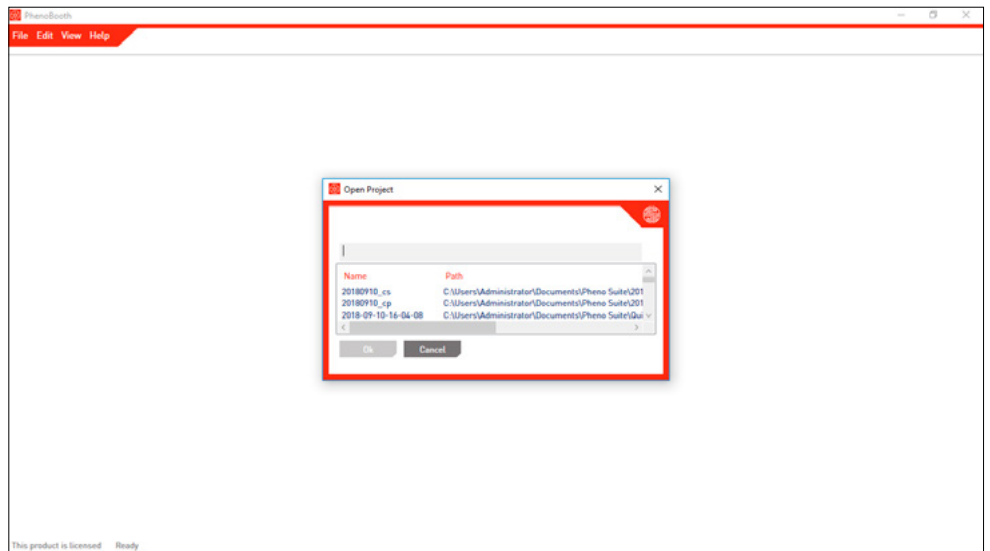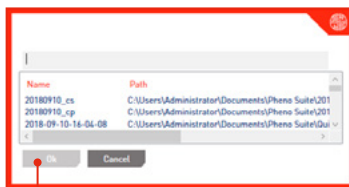

- If you want to open a previous project select Open Project and pick from the list of existing projects.
- Click OK to launch project.



COLONY SCREENING - IMAGE ACQUISITION

1

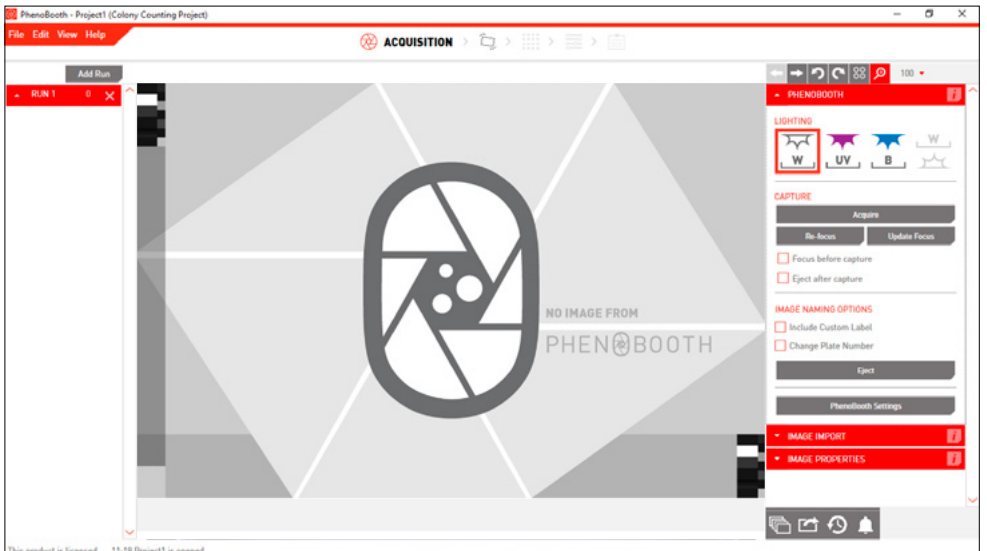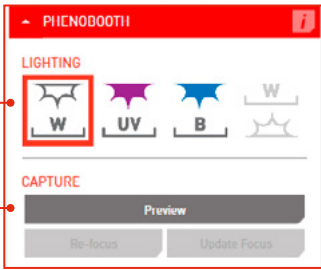- There are two options for acquiring an image: PhenoBooth+ or import an existing image.
- Select the desired method from the plug-in panel on the right of the screen. Press Preview to preview the image.



2

- If using the PhenoBooth+, click on that tab and select your desired light settings and select either Control or Experiment. Plate names are automatically generated.
- Insert a plate into PhenoBooth+ and press Acquire.
- Eject will allow you to remove the plate.
- Repeated until all of your plates have been captured.
- ➡ moves onto the pre-processing stage.

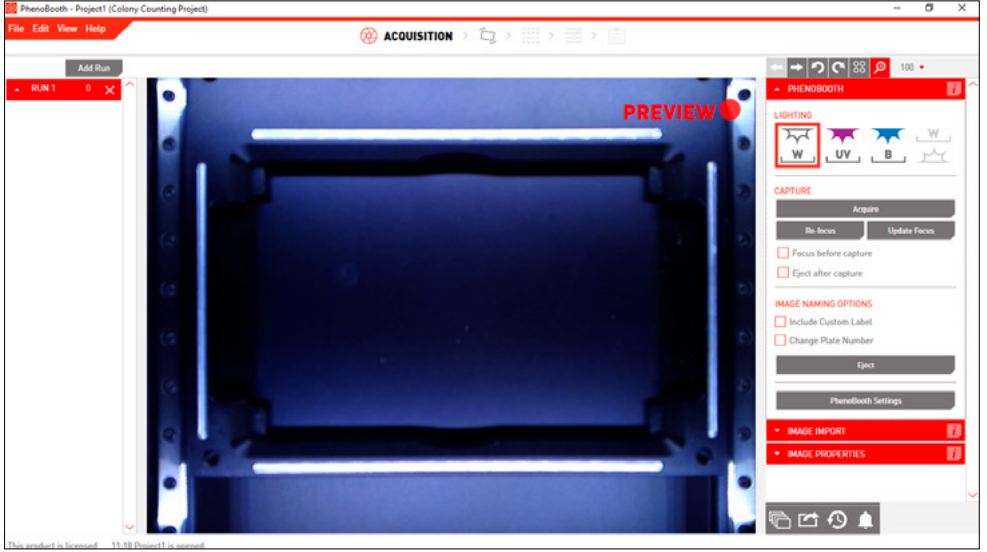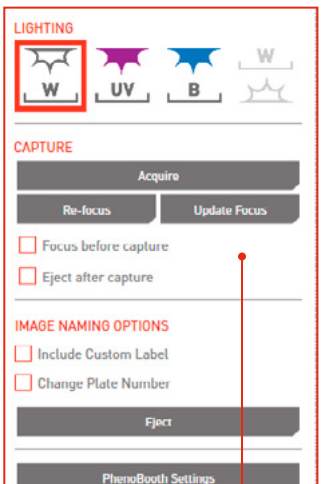
COLONY SCREENING - IMAGE ACQUISITION

3

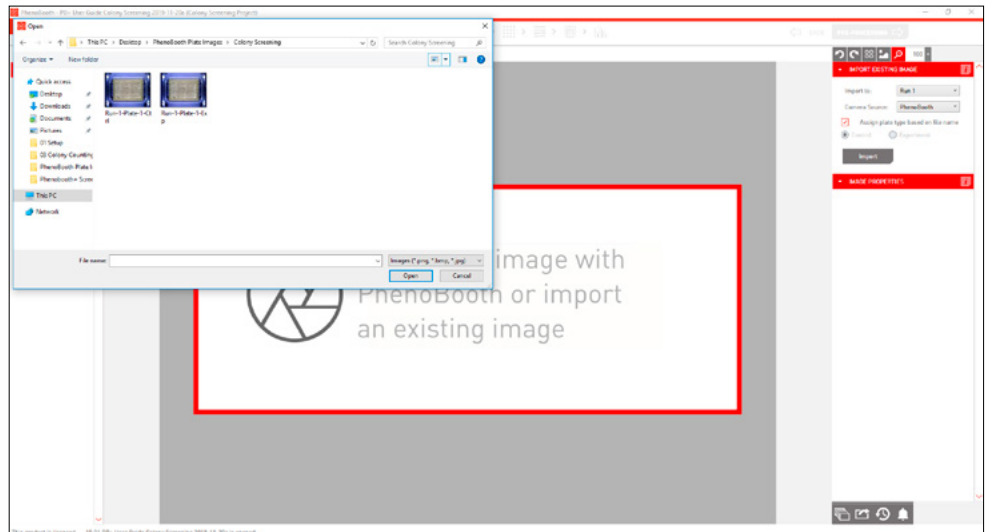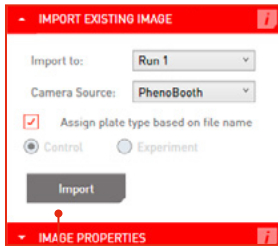

- If using existing images, click the Image Import tab. Here you can select which run the plate images will be associated. If this is your first run then Run 1 will be the only option. See p.49 to add new runs.
- Select whether the image is a control or experiment, or allow the PhenoBooth+ software to assign based on name. For example: *run-1-plate-1-ctrl* or *run-6-plate-9-exp*.
- Click Import to select images from your computer.


4

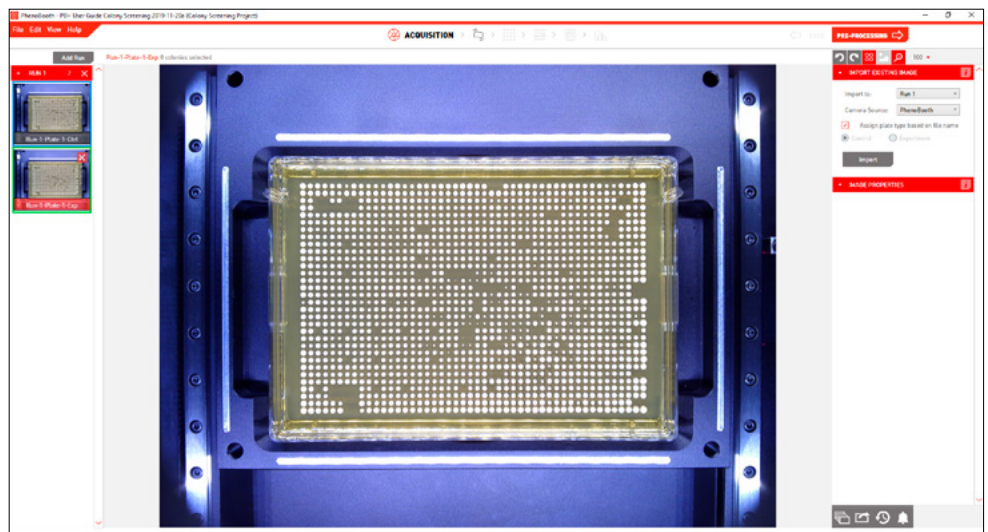

- Your captured or imported images now appear on the left of the screen. The currently selected plate is highlighted with red. Control plates have a blue border and experiment plates have a green border.
- Once all plates have been captured or imported, use the next stage button to move onto the pre-processing stage.



COLONY SCREENING - PRE-PROCESSING

5

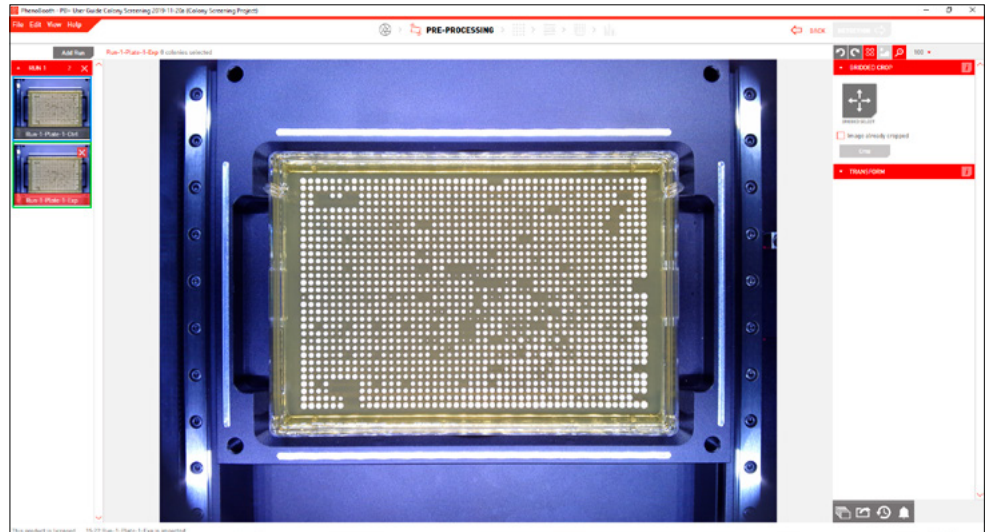

- Select which image to process first. If using the batch process option, this plate's settings will be applied to the rest of the plates in the run.




6

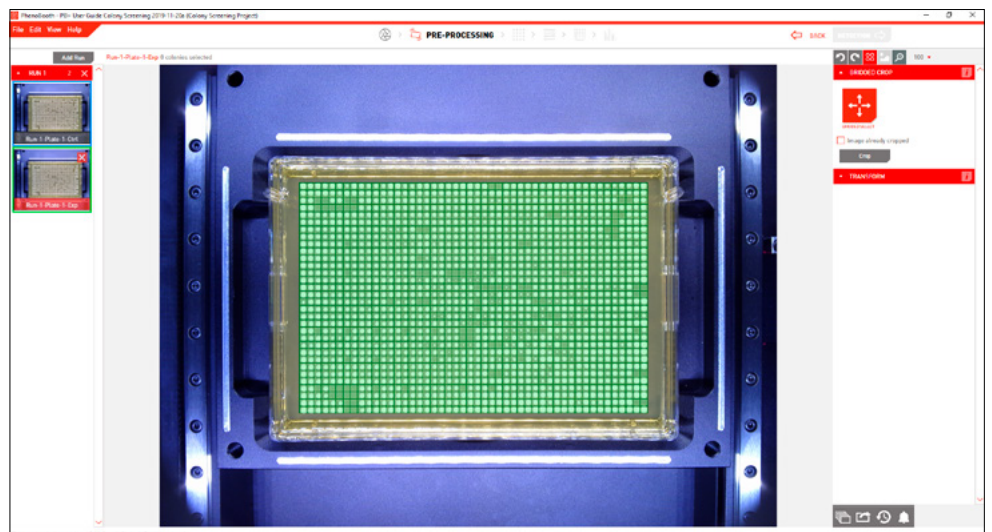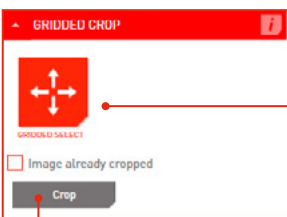

- In the Crop plug-in select Gridded Select.
- The grid will automatically overlay the plate image. You can also manually drag this grid over the plate with the mouse.
- Press Crop to remove the edge of the plate and beyond.
- Once cropped, you will automatically move to the detection stage.

COLONY SCREENING - DETECTION



7

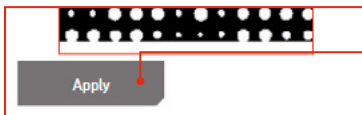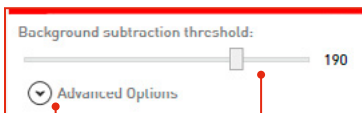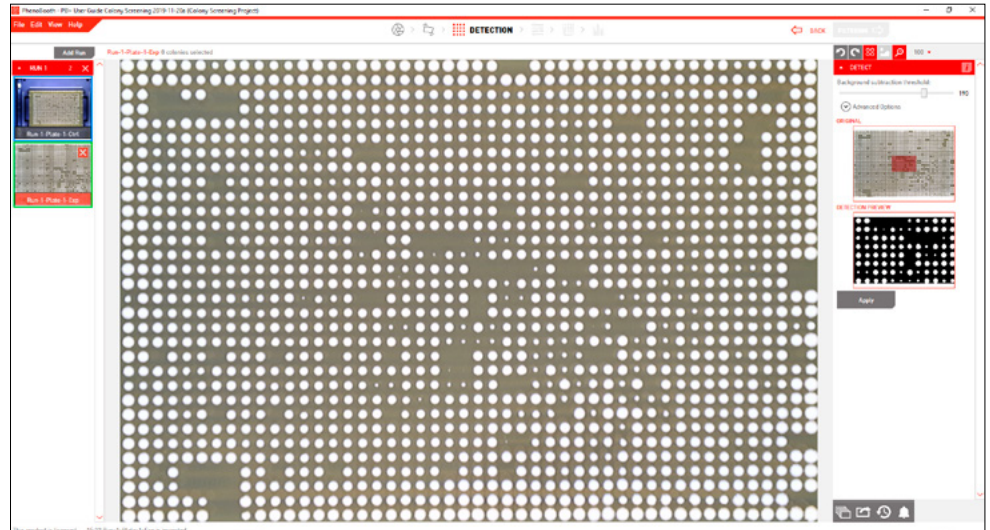


- In order to generate colony information, the background needs to be removed.
- Use the slider to adjust the filter intensity if necessary.
- If you have colonies that touch you can use the colony separation tool found in the advanced options.
- Check that all the smaller colonies are visible then click Apply.
- Once applied, you will automatically move to the filtering stage.


COLONY SCREENING - FILTERING

8

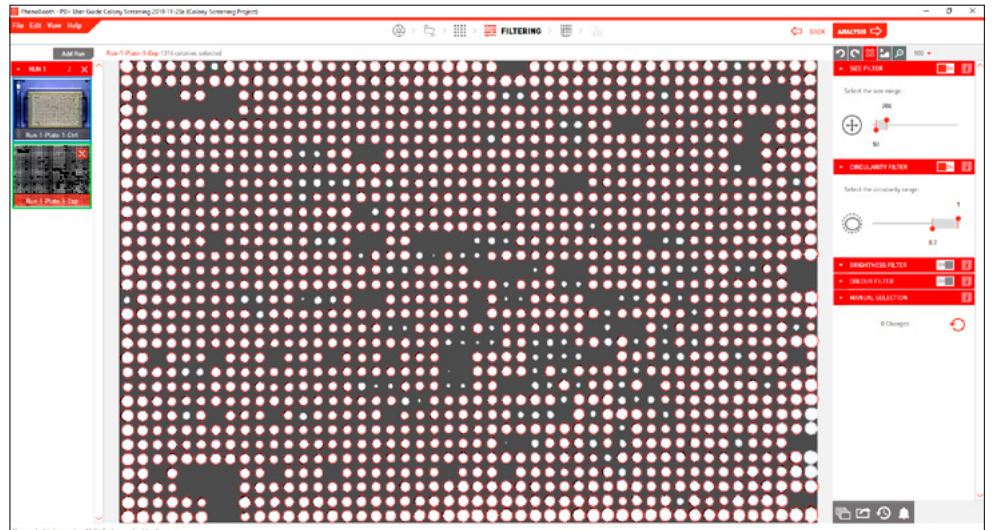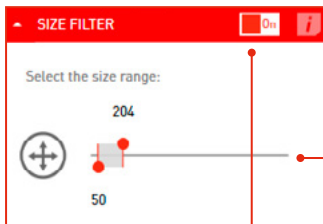

- Detected colonies are highlighted with a red circle.  turns highlights on or off.
- Use the filter options to set minimum and maximum values for size, circularity, brightness and colour for colony detection. You can use the size filter to discard items that are too large or small to be colonies.
- Turn filters on and off using the toggle button.
- Colonies can also be manually selected by clicking simply clicking on a colony to select or deselect it.


9

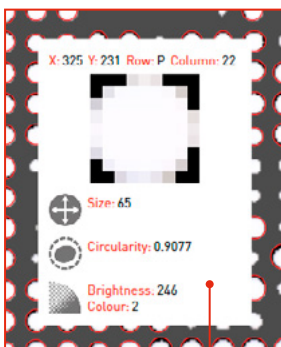

- Hover over a colony to see it's size, circularity and colour properties in the Colony Info plug-in tab.
- Now is the best time to batch process the rest of your plates with the same options using the instructions on p.47.
- Use the next stage button to move to the analysis stage.



COLONY SCREENING - ANALYSIS

10

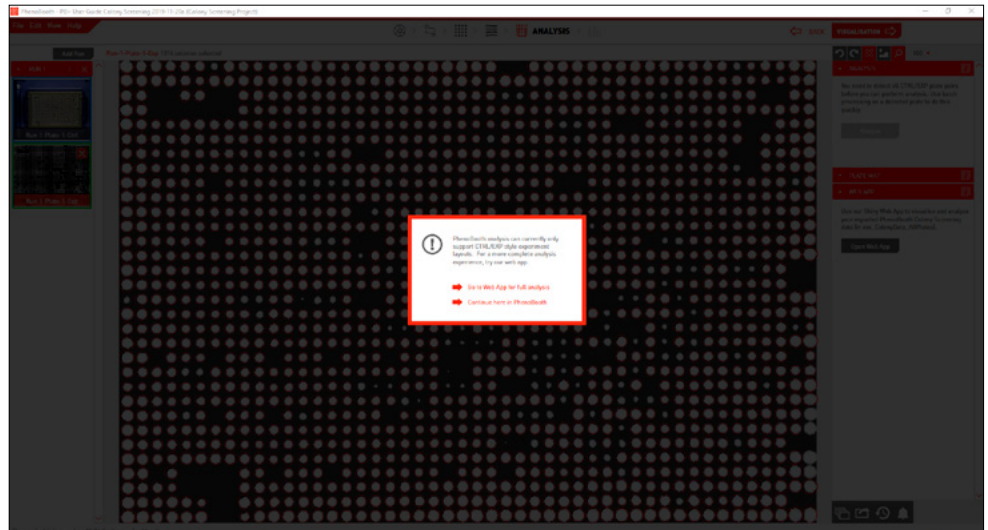

- The PhenoBooth+ software will handle simple control/experiment comparison. For more thorough analysis we have an online app.
- To move to the web app, select the Web App option from the pop-up window. You will be prompted to export your project at this stage.
- Click the Continue line to stay in PhenoBooth+.



11

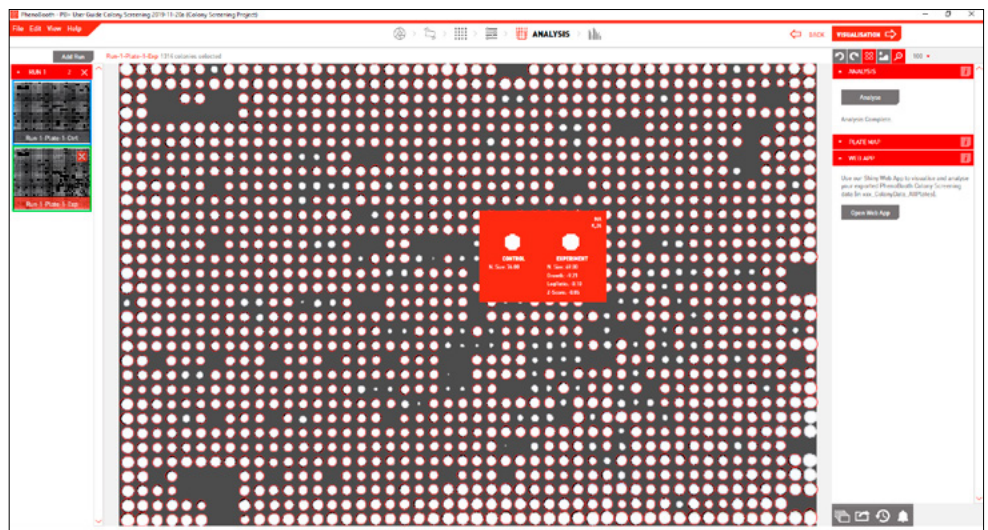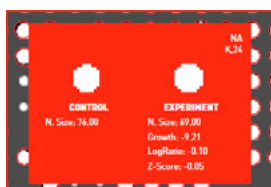

- To analyse colony growth difference between your current plate and it's control/experiment plate, select the Analysis tab and click Analyse.
- **NOTE:** To analyse you must first process all plates. To do this use the batch process instructions on p.47.
- Place the cursor over a colony to view the growth information for the control and experiment. Colony co-ordinates and gene information (if available) are shown in the top right corner of the info box. You can upload or select a plate map using the Plate Map plugin.
- Use the next stage button to move to the visualisation stage.

COLONY SCREENING - VISUALISATION

12

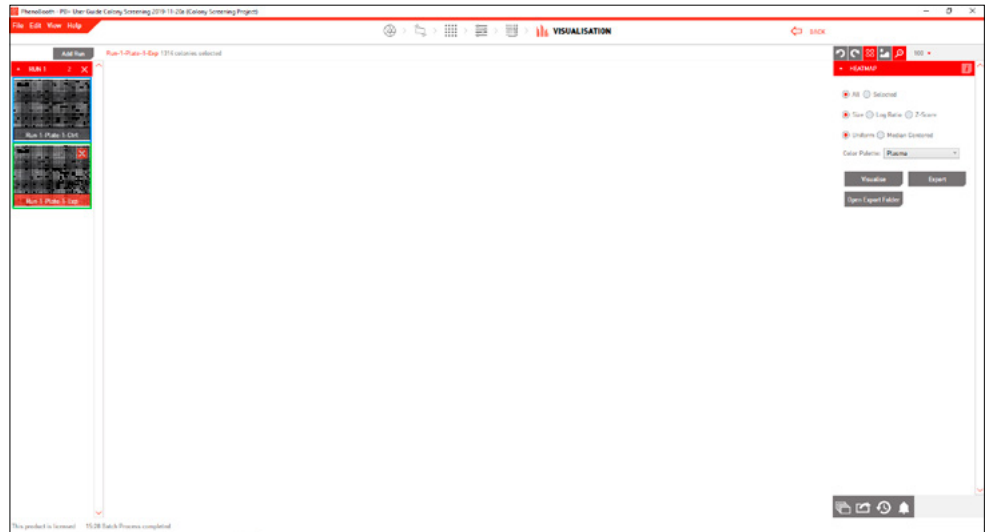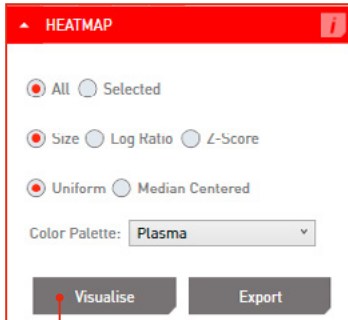

- To view heat maps of Size, Log Ratio or Z-score, choose your visualisation options from the Heat Map tab, then click Visualise.
- You can select to visualise all of your plates or just a selection.



13




- Boom! You can now see heat maps for your selected plates. Hover over the map to see the colony comparison info.

COLONY SCREENING - EXPORT

14

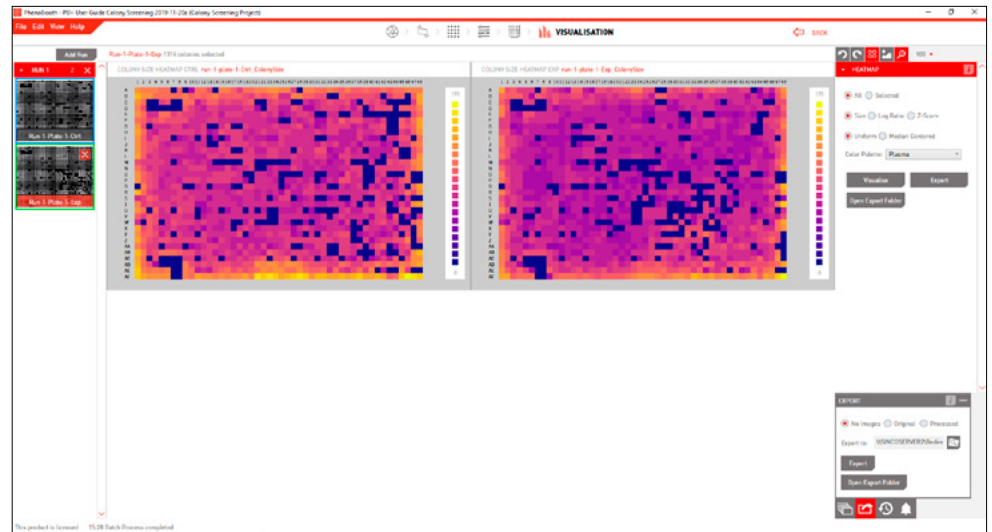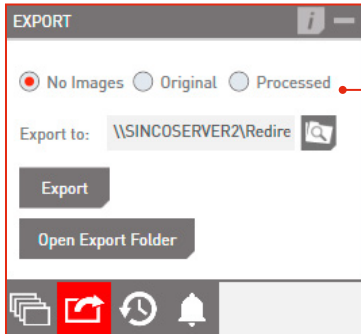

- The heatmap image can be exported in from the Heatmap plugin
- To export your entire project, use the Export tab. You can choose to export no images, processed images or the original images.
- Image formats available are: .bmp, .jpeg, .png or .tiff.
- The export includes: plate image (if selected), summary file, colony data table and colony data/analysis screen.
- The Summary File contains colony counts for all plates as well as colony imaging and filtering options information.
- The Colony Data Table contains position and properties of all detected colonies on all plates.

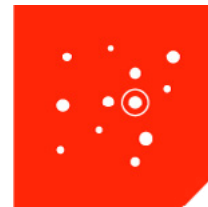


COLONY PICKING PROJECT

The Colony Picking package allows you to select individual colonies and export their coordinates to be picked using the Singer ROTOR. This section will guide you through the Colony Picking package for selection as well as the ROTOR software steps for picking.



There are 5 stages in the Colony Picking package: Image Acquisition, Pre-processing, Detection, Filtering and Stinger Export.

Start a new project then select Colony Picking.

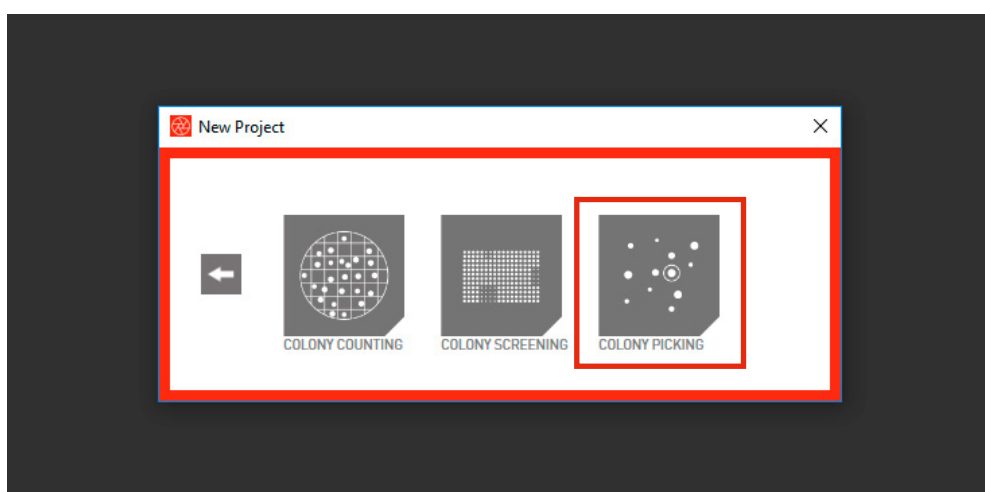



COLONY COUNTING - NEW PROJECT

1

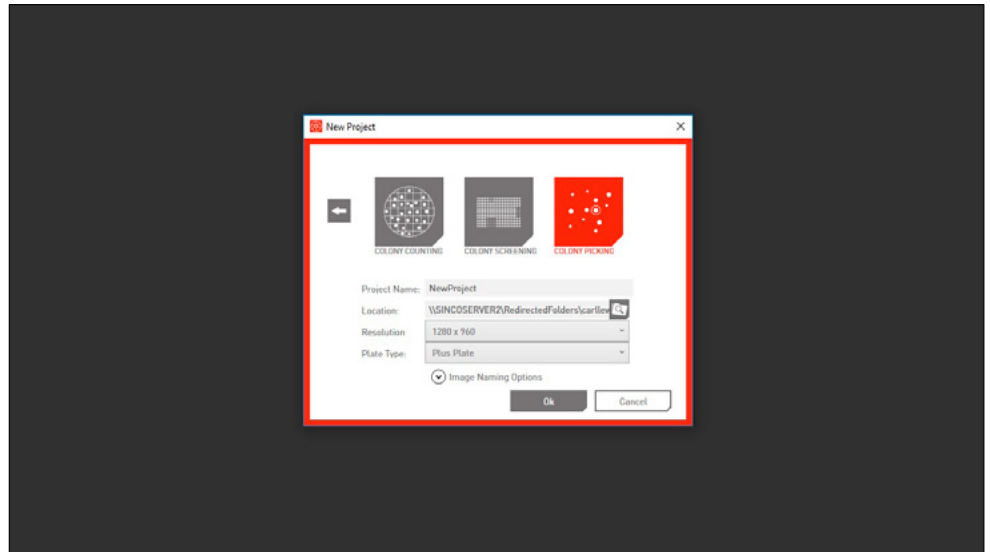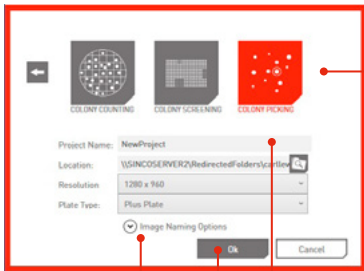

- Select the Colony Screening Icon. Icons will vary depending on which software packages you have purchased. Your chosen project type is highlighted in red.
- Name your project, select a save location and choose an image resolution and plate type.
- There are also a few optional naming settings you can use to customise your plate images.
- Click OK to launch project.



2

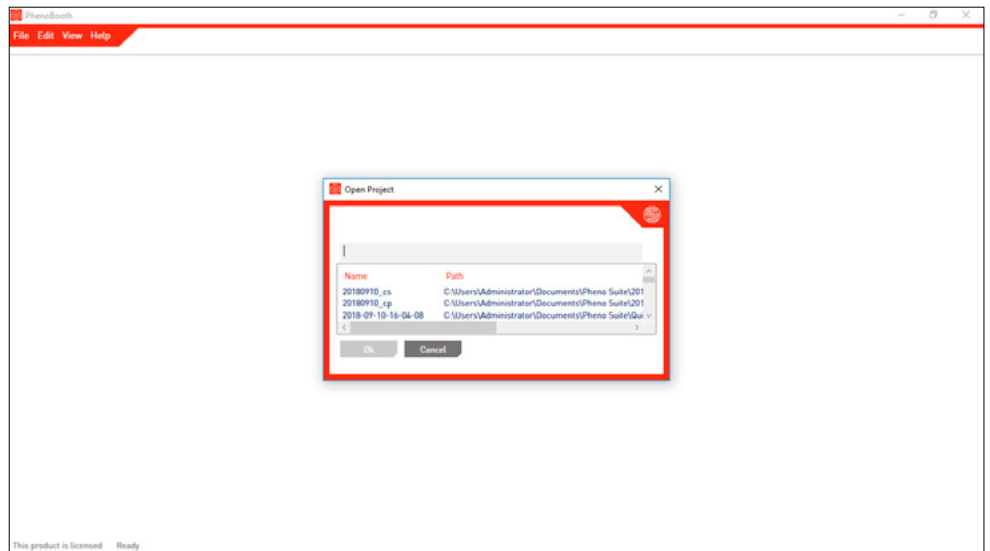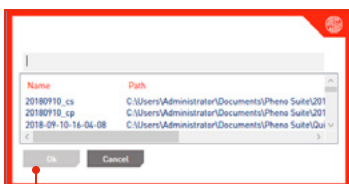

- If you want to open a previous project select Open Project and pick from the list of existing projects.
- Click OK to launch project.


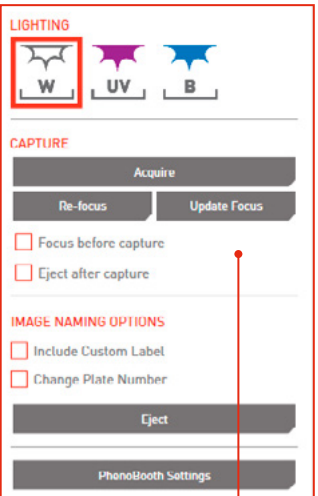
COLONY PICKING - IMAGE ACQUISITION

1

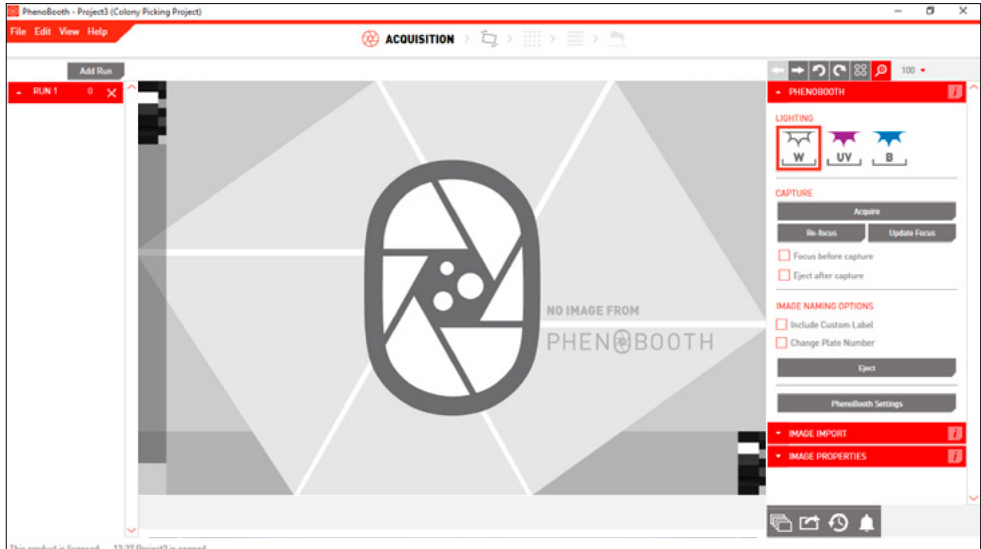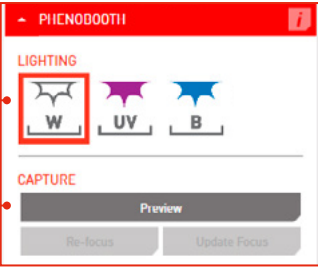- There are two options for acquiring an image: PhenoBooth+ or import an existing image. However, using imported images for colony picking is not recommended because the plate may have changed since the photo was taken, compromising accuracy.
- Select your lighting settings and press Preview to preview a live image from PhenoBooth+.


2

- If using the PhenoBooth+, click on that tab and select your desired light settings. Plate names are automatically generated.
- Insert a plate into PhenoBooth+ and press Acquire.
- Eject will allow you to remove the plate.
- Repeated until all of your plates have been captured.
- ➡ moves onto the pre-processing stage.

COLONY PICKING - IMAGE ACQUISITION



3

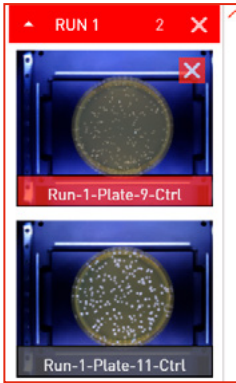

- Your captured images now appear on the left of the screen. The currently selected plate is highlighted with red.
- Use the next stage button to move to the pre-processing stage.


COLONY PICKING- PRE-PROCESSING & DETECTION

5

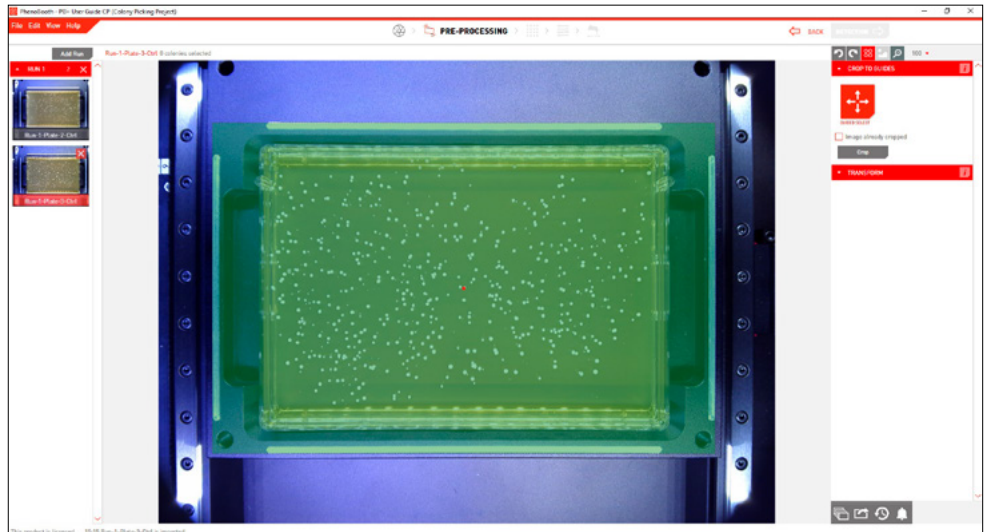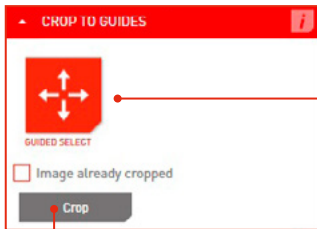

- In the Crop plug-in select Guided Select.
- The grid will automatically add a crop preview to the plate guides.
- Press Crop to remove unnecessary background.
- You will automatically move to the Detection stage.



6

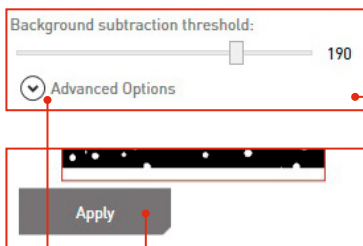


- The Background Subtraction plug-in is used to remove the plate background.
- Use the preview window to view colonies. Move the threshold bar if necessary to adjust the filter intensity.
- If you have colonies that touch you can use the colony separation tool found in the advanced options.
- When you're happy with the preview image, press Apply to remove the background.
- You will automatically move to the Analysis stage.


COLONY PICKING - FILTERING

7

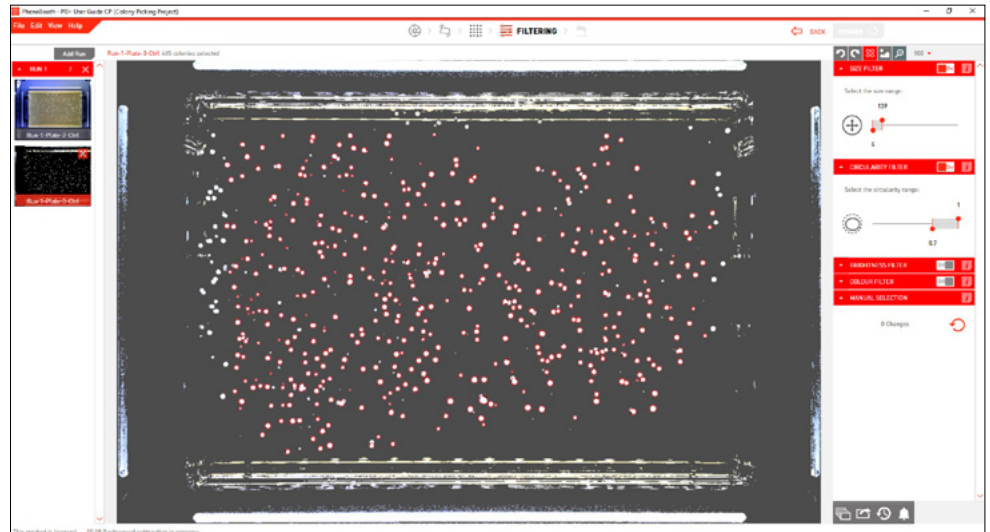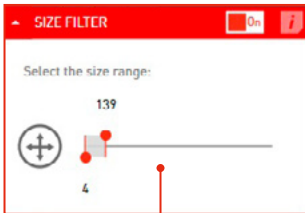

- Detected colonies are highlighted with a red circle.  turns highlights on or off.
- Use the filter options to set minimum and maximum values for size, circularity, brightness and colour for colony detection. You can use the size filter to discard items that are too large or small to be colonies.
- Colonies can also be manually selected by clicking on a colony to select or deselect it. Changes are displayed in the colony detection plugin tab.
- You can batch process the rest of your plates with the same options using the instructions on p.47.

8

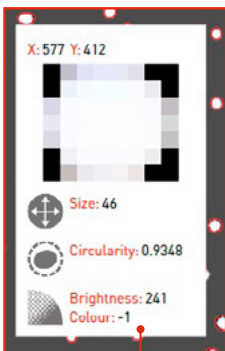


- Hover over a colony to see its size, circularity and colour properties in the Colony Info plug-in window.
- Once all of your plates have been processed (see batch process instructions on p.47) use the next stage button to move to the Stinger stage.


COLONY PICKING - STINGER EXPORT

9

Agar depth: 7

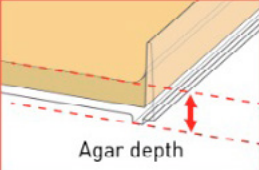


! Agar depth is the distance in millimetres between the bench and agar surface. This must be correct for accurate colony picking.

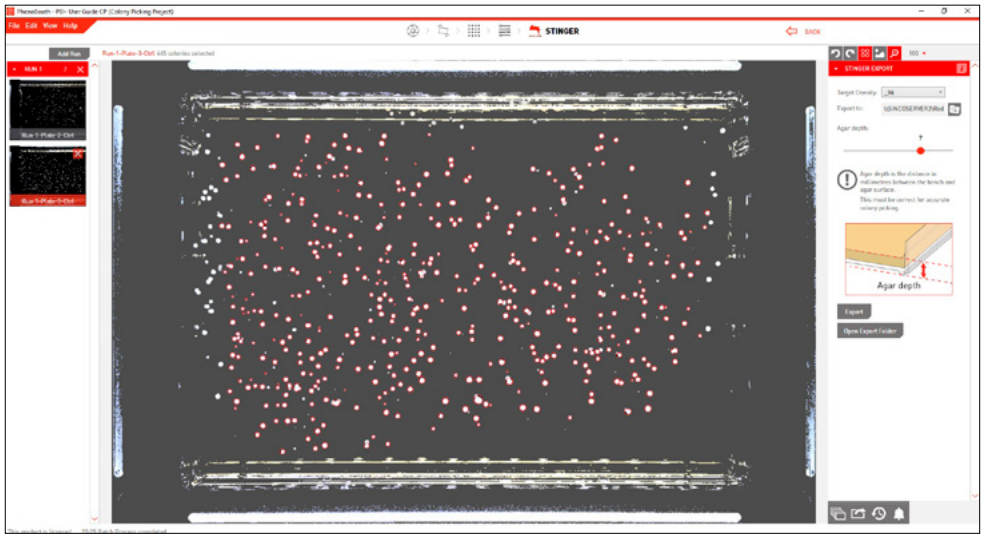


- Use the Stinger plugin to export colony locations for picking.
- Select the density you want the colonies arrayed in (96, 383, 1536 or 6144), select the file location.
- The depth of your agar will affect colony coordinates. Use the agar depth slider to set the correct depth.
- **NOTE:** Agar depth is the distance between the bench and the agar surface.
- The file will need to be saved to a USB stick so it can be transferred to the ROTOR HDA.

10

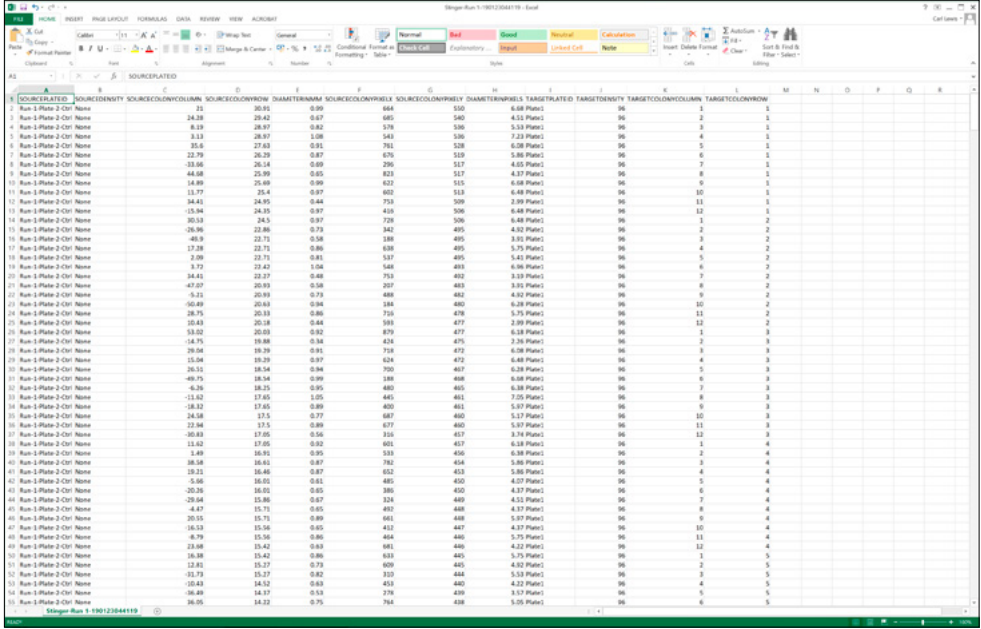

SOURCEID	SOURCEIDENITY	SOURCECOLONYNUMBER	SOURCECOLONYROW	DIAMETERMM	SOURCECOLONYRADIUS	SOURCECOLONYRADIUS	DIAMETERMM	TARGETIDENITY	TARGETCOLONYNUMBER	TARGETCOLONYROW
1	None	21	30.95	0.90	664	550	6.68 Plate1	96	2	1
2	Run 1-Plate 2-CH1-None	24.18	29.42	0.87	665	546	6.52 Plate1	96	2	1
3	Run 1-Plate 2-CH1-None	8.13	28.97	0.82	578	536	5.53 Plate1	96	3	1
4	Run 1-Plate 2-CH1-None	3.13	28.97	1.08	543	536	7.22 Plate1	96	4	1
5	Run 1-Plate 2-CH1-None	35.6	27.63	0.91	761	528	6.58 Plate1	96	5	1
6	Run 1-Plate 2-CH1-None	22.79	26.29	0.87	676	519	5.86 Plate1	96	6	1
7	Run 1-Plate 2-CH1-None	13.66	26.14	0.95	276	517	4.65 Plate1	96	7	1
8	Run 1-Plate 2-CH1-None	44.68	25.99	0.85	823	517	4.37 Plate1	96	8	1
9	Run 1-Plate 2-CH1-None	14.48	25.48	0.99	612	515	6.48 Plate1	96	9	1
10	Run 1-Plate 2-CH1-None	13.77	25.4	0.97	602	513	6.48 Plate1	96	10	1
11	Run 1-Plate 2-CH1-None	34.41	24.95	0.84	734	509	2.99 Plate1	96	11	1
12	Run 1-Plate 2-CH1-None	15.94	24.95	0.97	419	506	6.48 Plate1	96	12	1
13	Run 1-Plate 2-CH1-None	30.13	24.5	0.97	726	506	6.48 Plate1	96	1	2
14	Run 1-Plate 2-CH1-None	28.96	23.86	0.73	842	495	4.92 Plate1	96	2	2
15	Run 1-Plate 2-CH1-None	48.9	22.71	0.58	338	495	3.91 Plate1	96	3	2
16	Run 1-Plate 2-CH1-None	17.28	22.71	0.86	638	495	5.75 Plate1	96	4	2
17	Run 1-Plate 2-CH1-None	2.89	22.71	0.81	537	495	5.41 Plate1	96	5	2
18	Run 1-Plate 2-CH1-None	3.77	22.42	1.04	541	483	6.96 Plate1	96	6	2
19	Run 1-Plate 2-CH1-None	34.41	22.27	0.48	753	482	5.13 Plate1	96	7	2
20	Run 1-Plate 2-CH1-None	47.07	20.93	0.58	207	483	3.91 Plate1	96	8	2
21	Run 1-Plate 2-CH1-None	5.11	20.93	0.73	688	482	4.92 Plate1	96	9	2
22	Run 1-Plate 2-CH1-None	56.49	20.63	0.94	344	480	6.28 Plate1	96	10	2
23	Run 1-Plate 2-CH1-None	28.75	20.33	0.88	716	478	5.75 Plate1	96	11	2
24	Run 1-Plate 2-CH1-None	10.43	20.18	0.44	589	477	2.99 Plate1	96	12	2
25	Run 1-Plate 2-CH1-None	53.02	20.03	0.92	879	477	6.18 Plate1	96	1	3
26	Run 1-Plate 2-CH1-None	14.75	19.88	0.84	414	475	2.76 Plate1	96	2	3
27	Run 1-Plate 2-CH1-None	29.04	19.29	0.91	718	472	6.28 Plate1	96	3	3
28	Run 1-Plate 2-CH1-None	15.04	19.29	0.97	614	472	6.48 Plate1	96	4	3
29	Run 1-Plate 2-CH1-None	26.11	18.54	0.94	700	467	6.28 Plate1	96	5	3
30	Run 1-Plate 2-CH1-None	48.75	18.54	0.99	339	468	6.68 Plate1	96	6	3
31	Run 1-Plate 2-CH1-None	4.26	18.25	0.95	480	465	6.18 Plate1	96	7	3
32	Run 1-Plate 2-CH1-None	13.62	17.85	1.05	465	461	7.05 Plate1	96	8	3
33	Run 1-Plate 2-CH1-None	22.82	17.85	0.89	609	461	5.97 Plate1	96	9	3
34	Run 1-Plate 2-CH1-None	24.58	17.5	0.77	687	460	5.17 Plate1	96	10	3
35	Run 1-Plate 2-CH1-None	22.94	17.5	0.89	617	460	5.97 Plate1	96	11	3
36	Run 1-Plate 2-CH1-None	36.83	17.05	0.56	316	457	3.74 Plate1	96	12	3
37	Run 1-Plate 2-CH1-None	13.62	17.05	0.92	601	457	6.18 Plate1	96	1	4
38	Run 1-Plate 2-CH1-None	1.49	16.91	0.95	519	456	6.18 Plate1	96	2	4
39	Run 1-Plate 2-CH1-None	38.38	16.43	0.87	782	454	5.86 Plate1	96	3	4
40	Run 1-Plate 2-CH1-None	19.11	16.46	0.87	613	453	5.86 Plate1	96	4	4
41	Run 1-Plate 2-CH1-None	5.46	16.02	0.81	485	450	4.97 Plate1	96	5	4
42	Run 1-Plate 2-CH1-None	20.26	16.02	0.85	386	450	4.97 Plate1	96	6	4
43	Run 1-Plate 2-CH1-None	29.04	15.86	0.87	324	449	4.51 Plate1	96	7	4
44	Run 1-Plate 2-CH1-None	4.67	15.71	0.85	492	448	4.97 Plate1	96	8	4
45	Run 1-Plate 2-CH1-None	30.15	15.71	0.89	661	448	5.97 Plate1	96	9	4
46	Run 1-Plate 2-CH1-None	16.13	15.56	0.85	412	447	4.37 Plate1	96	10	4
47	Run 1-Plate 2-CH1-None	6.79	15.56	0.86	464	446	5.75 Plate1	96	11	4
48	Run 1-Plate 2-CH1-None	23.88	15.42	0.83	681	446	4.22 Plate1	96	12	4
49	Run 1-Plate 2-CH1-None	38.38	15.42	0.86	633	445	5.75 Plate1	96	1	5
50	Run 1-Plate 2-CH1-None	12.81	15.27	0.79	609	445	4.82 Plate1	96	2	5
51	Run 1-Plate 2-CH1-None	33.73	15.27	0.82	313	444	5.53 Plate1	96	3	5
52	Run 1-Plate 2-CH1-None	10.43	14.12	0.83	613	440	4.52 Plate1	96	4	5
53	Run 1-Plate 2-CH1-None	36.49	14.17	0.53	278	439	3.17 Plate1	96	5	5
54	Run 1-Plate 2-CH1-None	36.05	14.12	0.79	764	438	5.05 Plate1	96	6	5

- The exported .CSV file contains all information on the selected colonies.
- It's now time to load this file into the ROTOR HDA.

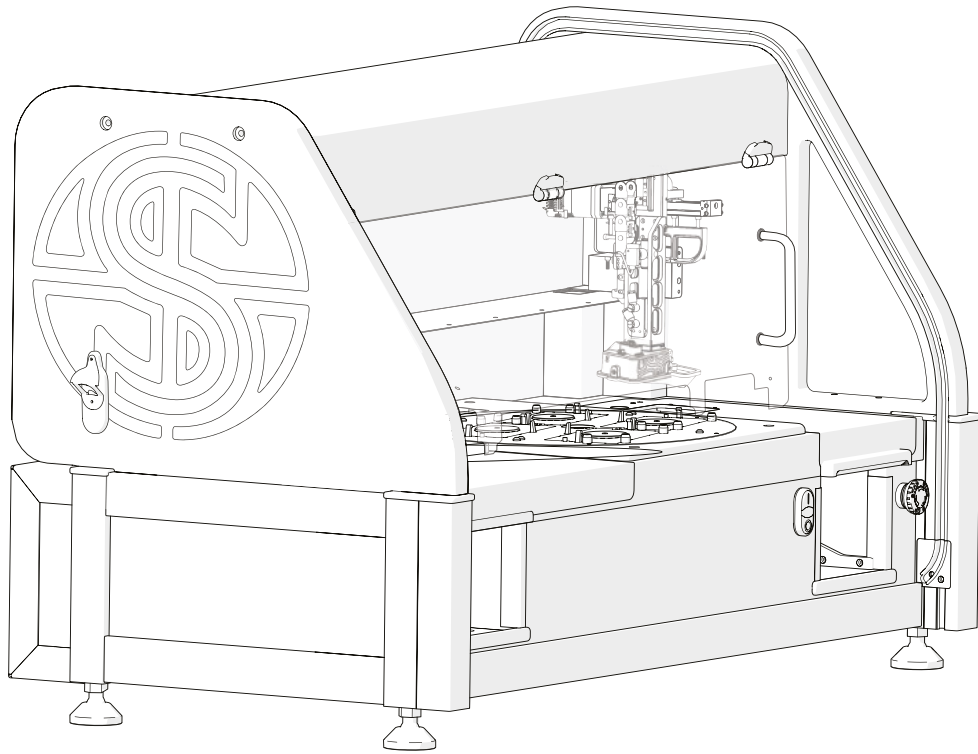


Now head over to the ROTOR HDA.

Please make sure The Stinger single colony picker is installed.


1

- Turn on the ROTOR and ensure that The Stinger is installed.
- On the ROTOR start-up screen you will be asked to select an operation mode. Select the Programming /Stinger option.

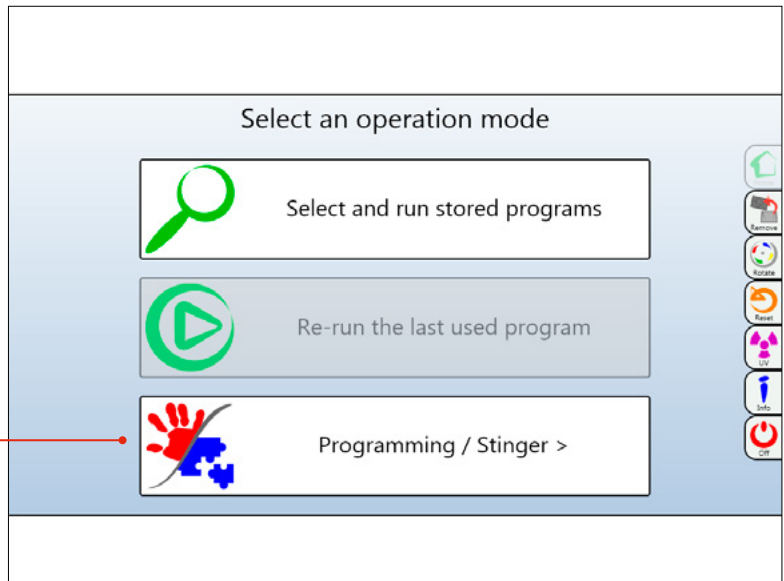


2

- Select the option labelled Stinger.

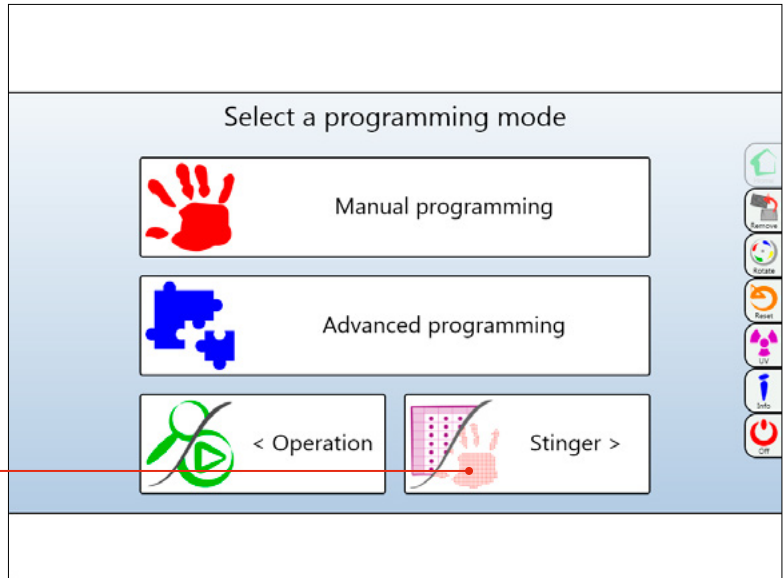

3


- On the Stinger Home Screen, select the Import Spreadsheet option.

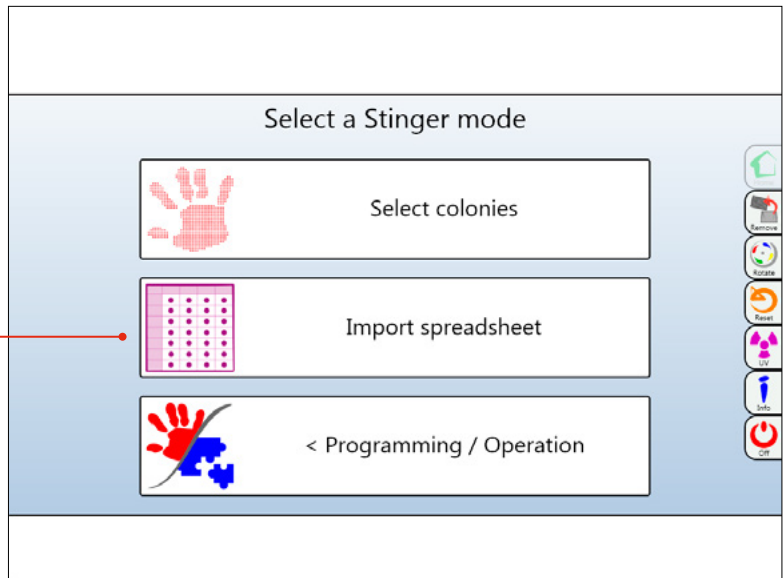

ROTOR HDA / THE STINGER

4

- This screen will appear if no USB memory stick is inserted. Insert your USB memory stick now.

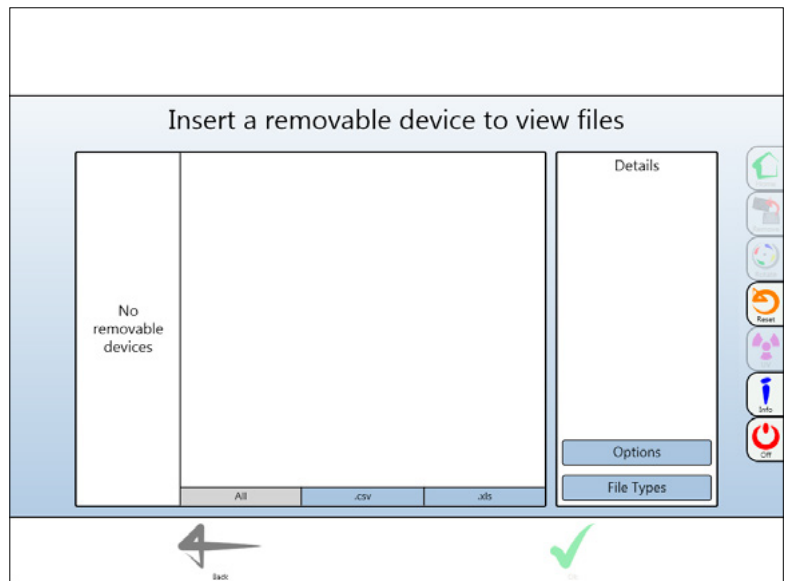


5

- Once your USB memory stick has been inserted, all .csv files in the root directory are displayed.

Memory stick

Files available

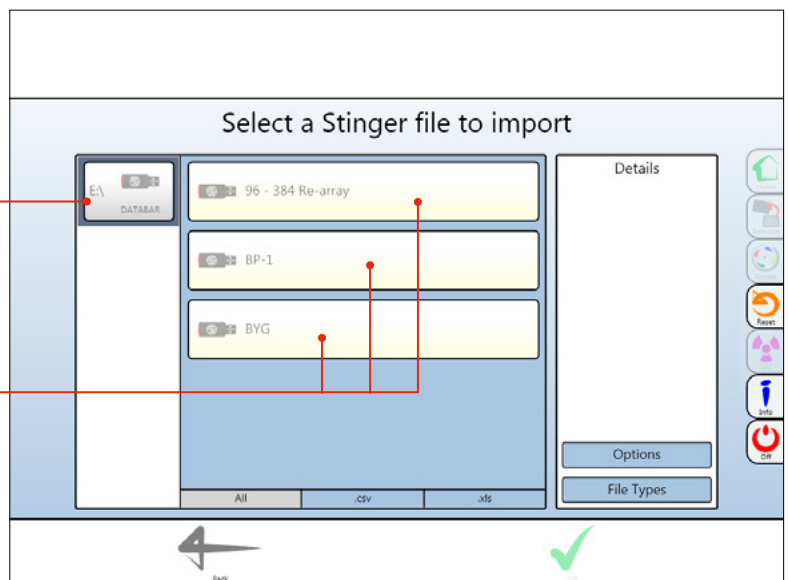


6

- Select the file you exported from PhenoSuite.

- Details of the file are displayed on the right.


- Select OK to run chosen file.

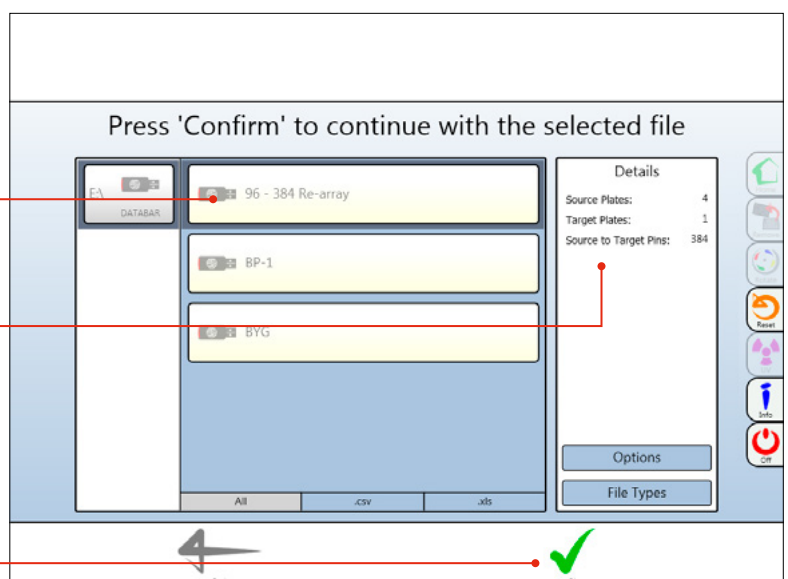

ROTOR HDA / THE STINGER

7


- The final confirmation screen will appear. Click on the tabs to change advanced options if necessary.

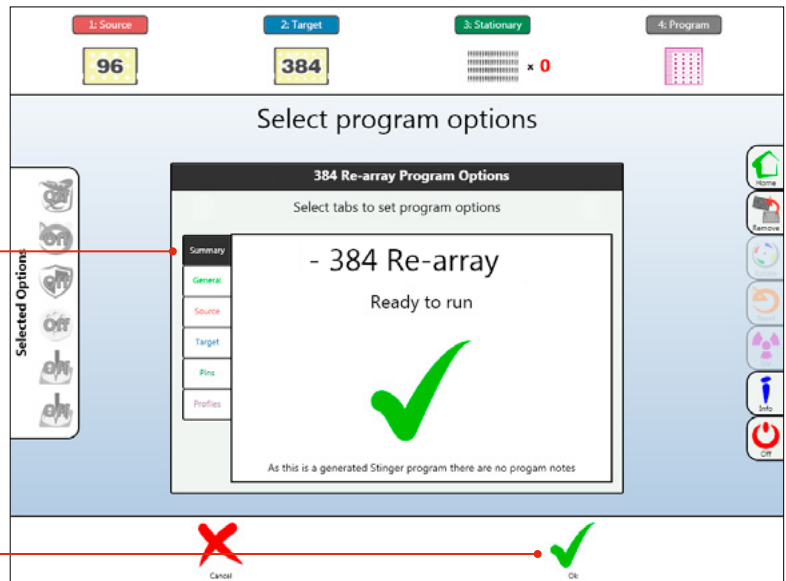


- Select OK to start program.

8


- Live program information will be updated throughout the operation.
- Program: resume, pause, stop / abort.

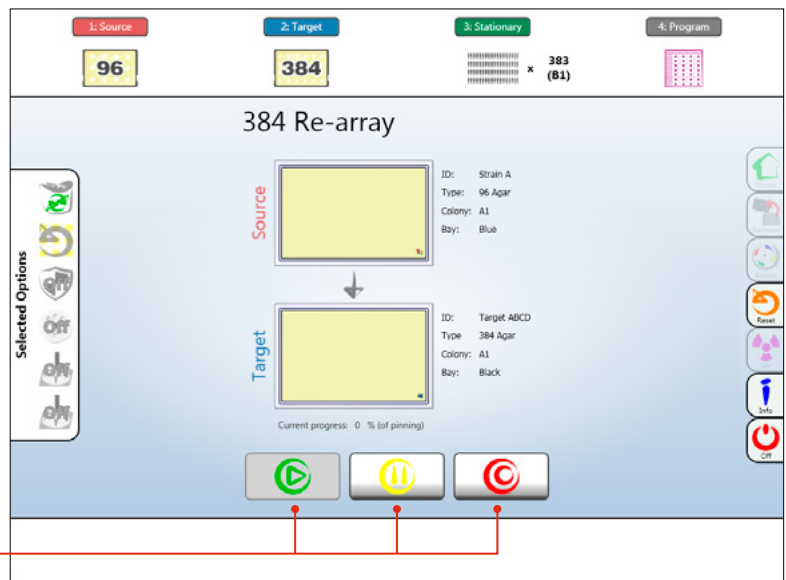

SOURCE PLATE

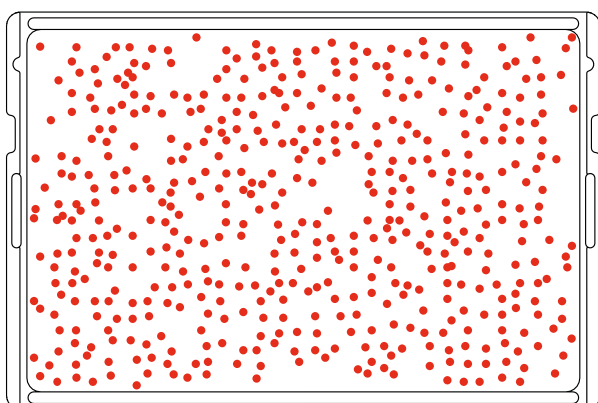

TARGET PLATE

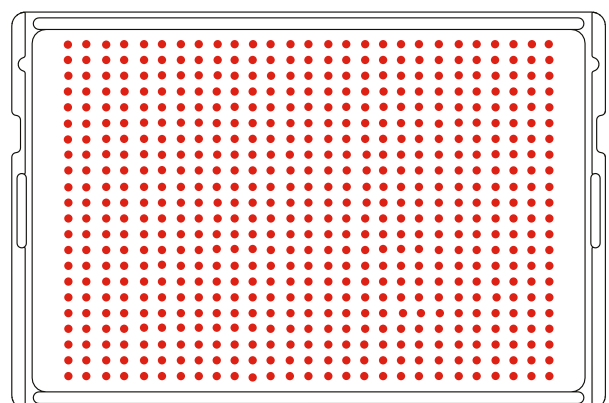

BEST PRACTICE

We want you to have a trouble-free time with your PhenoBooth+ so we've assembled some handy hints, tips and troubleshooting answers to keep you on the straight and narrow.

SOFTWARE ICONS

GENERAL TOOLS



UNDO. Undo the last process completed.




REDO. If you just used the undo tool, you can redo that process again with this button.



HIGHLIGHT COLONIES. Turn on or off detected colony highlights.



ZOOM. Use to zoom in on colonies of interest.

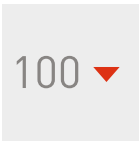

ZOOM PERCENTAGE. Selects zoom percentage of current plate.


STAGE NAVIGATION



IMAGE ACQUISITION.



ANALYSIS. Colony Screening package only.

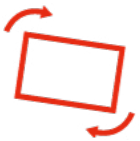

PRE-PROCESSING.



VISUALISATION. Colony Screening package only.



DETECTION.



STINGER. Colony Picking package only.



FILTERING.

SOFTWARE ICONS

NOTIFICATIONS



CONNECT PHENOBOOTH+. PhenoBooth+ needs to be connected.



CONNECTION PROBLEM. There is a problem with the PhenoBooth+ connection. Try reconnecting the USB 3.0 cable less slowly, see p.2.



LIGHTING TIMED OUT. If left for too long the lighting goes on standby.



TRAY CRASH. The input tray has crashed. Ensure plate is loaded correctly and the lid has been removed.



CAPTURING. PhenoBooth+ is taking an image.



TRAY ACTIVE. Input tray is moving.

LIGHTING OPTIONS



WHITE TOP LIGHT. White, flat-field incident lighting.



UV TOP LIGHT. Ultra-violet, flat-field incident lighting.




BLUE TOP LIGHT. Blue, flat-field incident lighting.




WHITE BOTTOM LIGHT. White, flat-field, transmitted light.

FEEDBACK LIGHT



 GREEN. All okay

 FLASHING RED FAST. Tray crash, Please wait while the tray automatically re-homes.

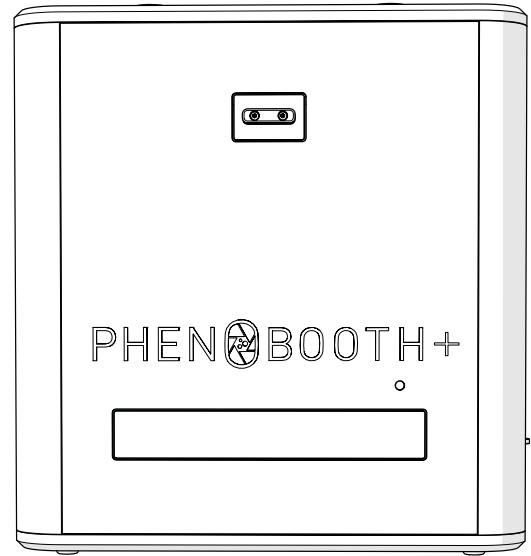


BATCH PROCESSING

1

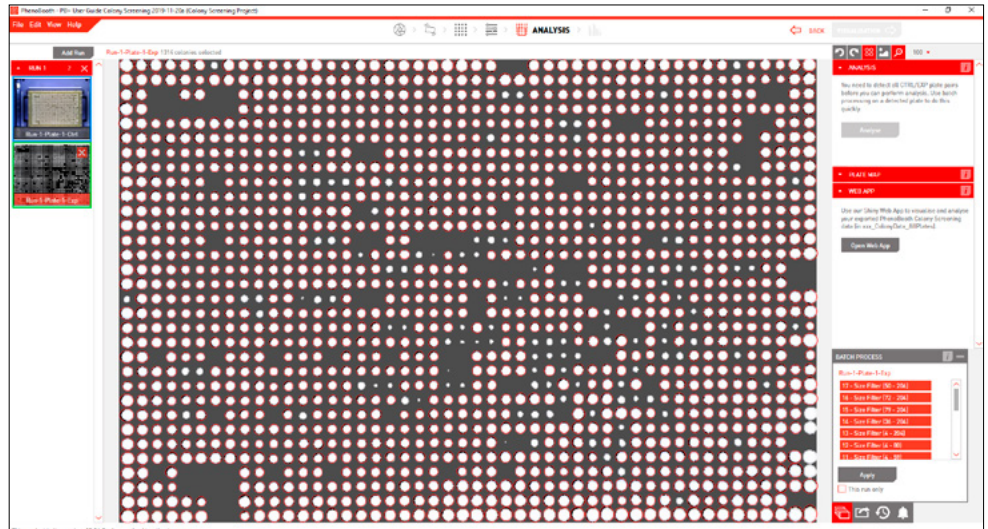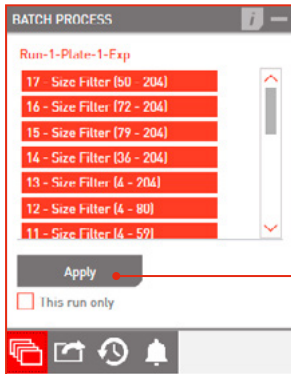

- Using the same settings for all plate images in a run will not only increase the speed of analysis but also ensure consistent and reliable information. A good time to batch process all of your plates is after all filtering has taken place.
- Click the Batch Process tab all processes are selected by default. Deselect any process you want to exclude.
- Click Apply and batch processing will begin. This may take a while for larger numbers of images.



2

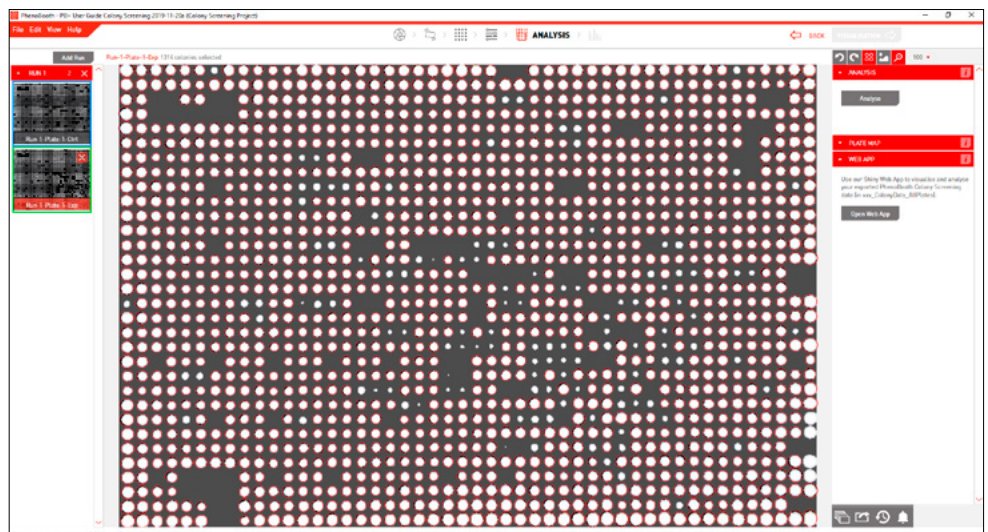


- All plate images in your current run have now been processed with the same options as your original image.
- You can see previews of the rest of your plates on the left of the screen. Click a plate to see a full screen view.


HISTORY PLUG-IN

1

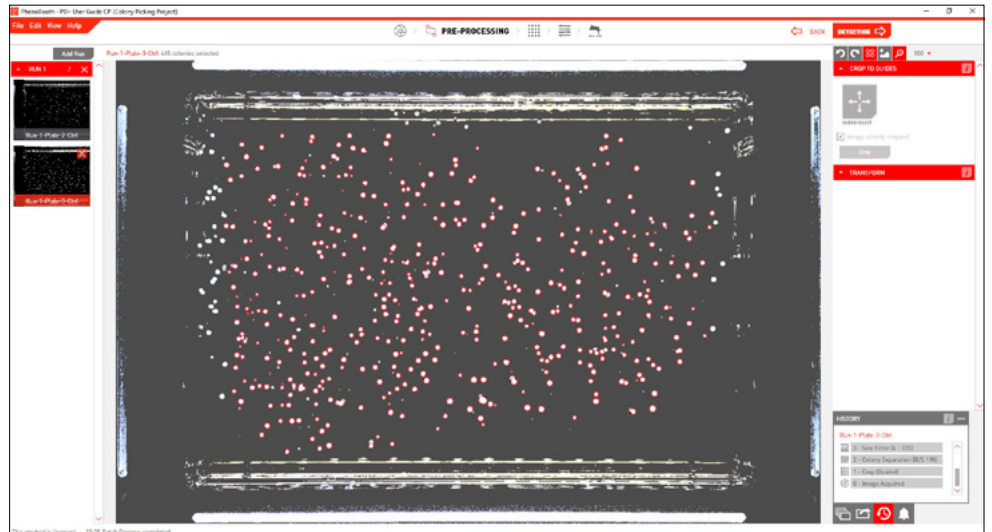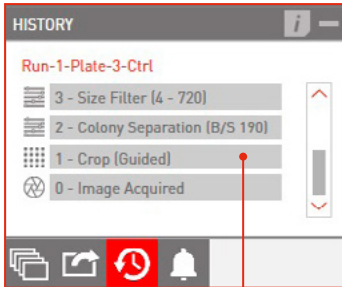

- If you want to undo the last change you made, simply click the undo button: 
- If you want to go back further then use the History plug-in. Open the History plug-in tab and select the process you want to revert back to. The list is in chronological order with the most recent process at the top.
- In this example we're reverting back to when we cropped the image.

2

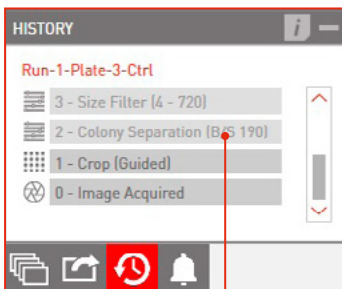

- Now, the last process applied to the image is crop. All processes after crop have been discarded.
- Discarded processes are marked as light grey in the History plug-in. You can revisit any of these processes by clicking on them, until you make a change to the plate image outside of the History plug-in.


ADD NEW PLATE RUN

1



- Your plates can be organised into separate runs for different groups of experiments.
- To add a new run, simply click Add Run in the Plate Preview section on the left of the screen.


2

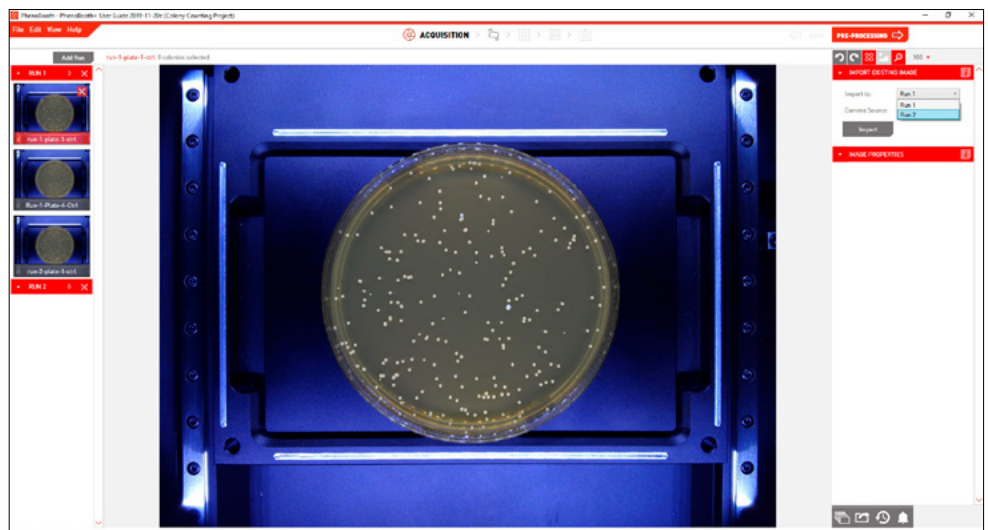

- When acquiring images, whether capturing using a PhenoBooth+ or importing existing images, you'll be asked to select which run you want your images to be filed under.
- Here you will see all the runs you have added so far.



IMPORT / EXPORT PROJECTS

1

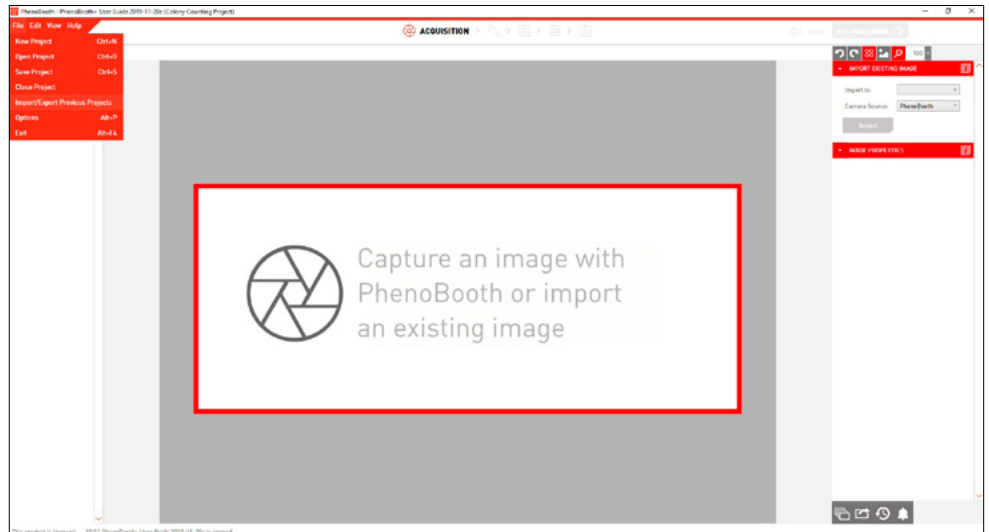

- PhenoBooth+ allows you to move projects from computer to computer. To do this, select the Import/Export Projects option in the File menu.


2

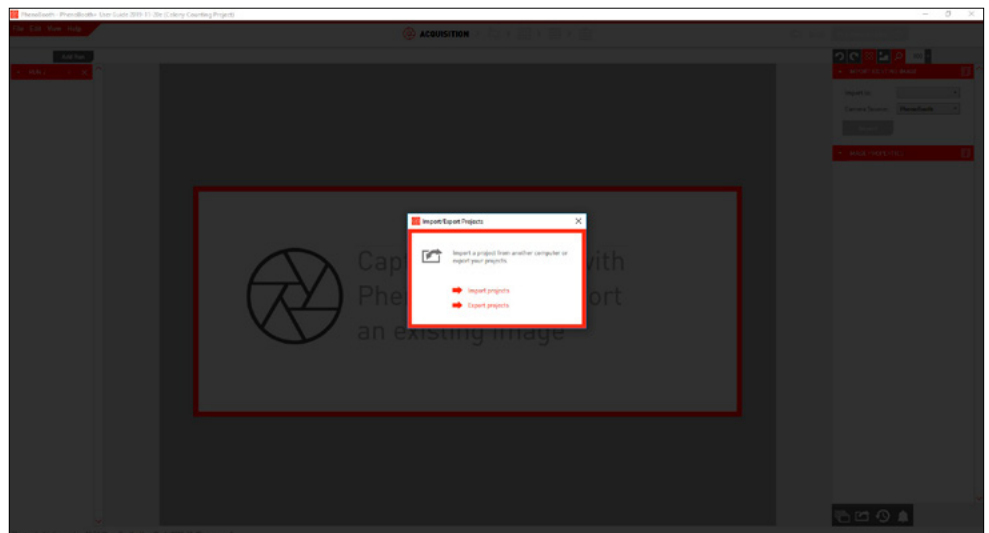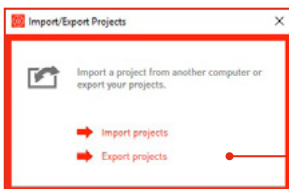

- In the pop-up window, select whether you want to import or export projects.


IMPORT / EXPORT PROJECTS

3

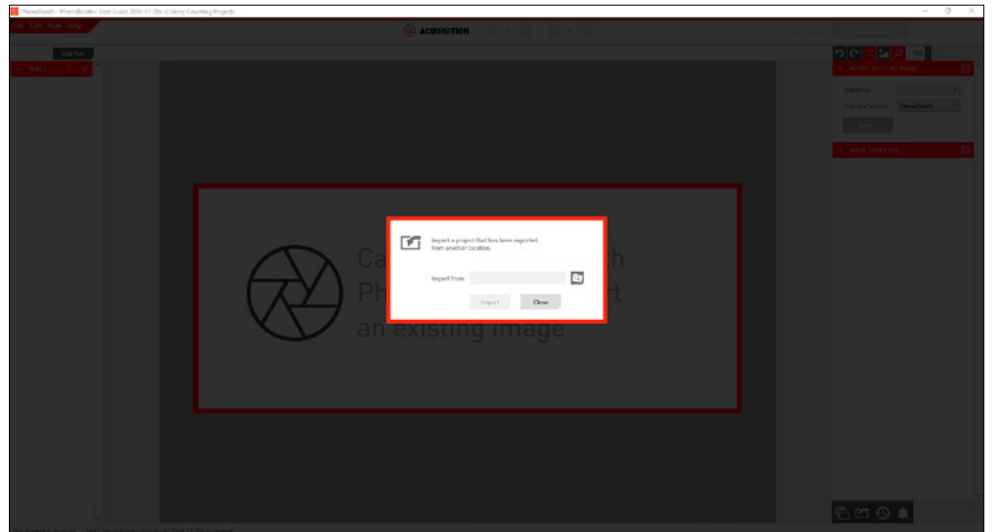


- For exporting, select the export location and then press Export.

4

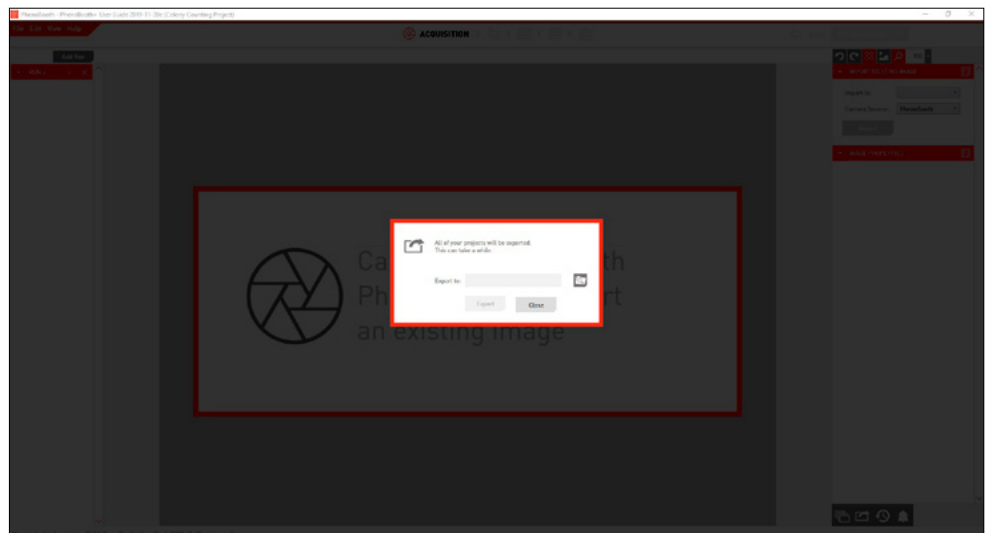


- For importing, select the import location and then press Import.


USING FILTERS

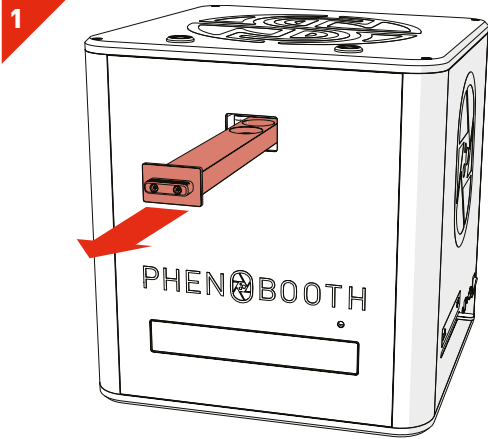


- Remove filter tray fully.

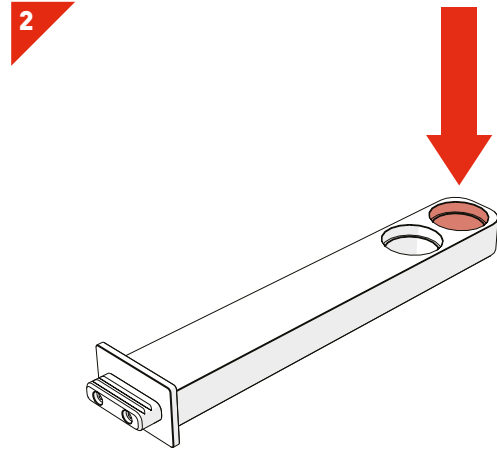

- Place the desired filter into the end filter slot.
- The PhenoBooth+ accepts standard 25mm circular filters with a maximum height of 5mm.

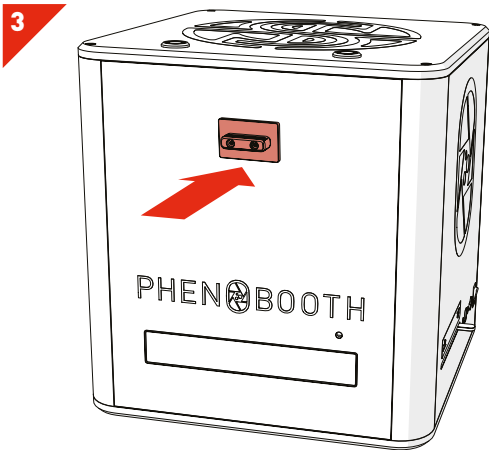


- Reinsert the filter tray fully until you feel the click. This will set the tray to the no filter position.

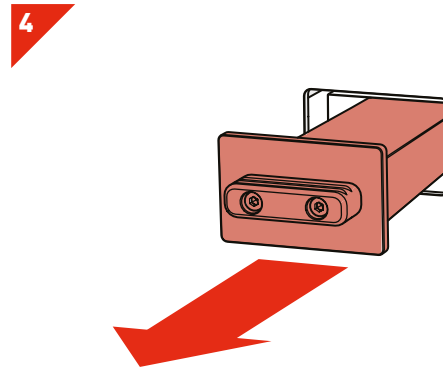


- To select the filter for use, pull the tray out slightly to the filter position — you will feel it click into place.

SETTING AGAR HEIGHT

1

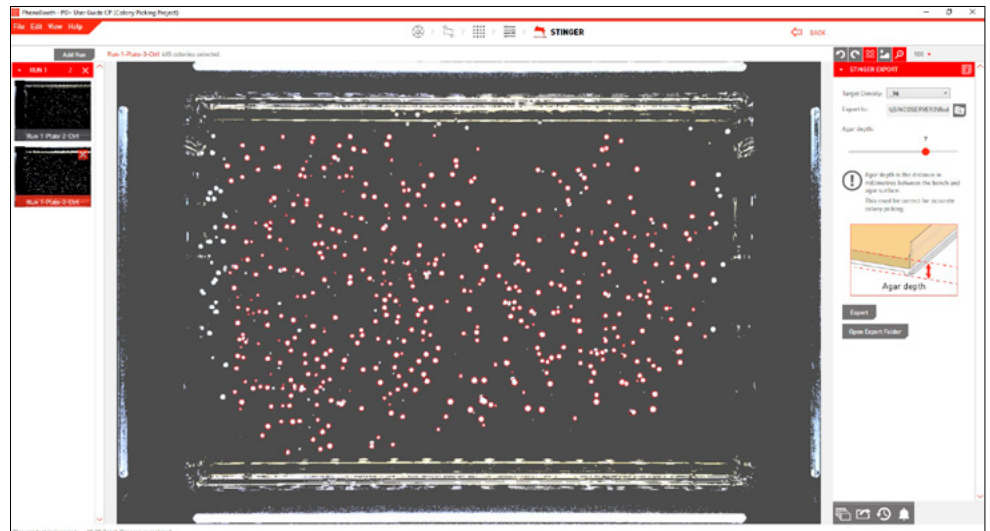

- The PhenoBooth+ is capable of using a variety of different plates which can have different base thicknesses. To account for this, we ask users to set their agar height before exporting to The Stinger colony picker.
- Agar height is measured as the distance from the top of the agar (trying to ignore the meniscus) to the bottom of the plate. See the diagram below.

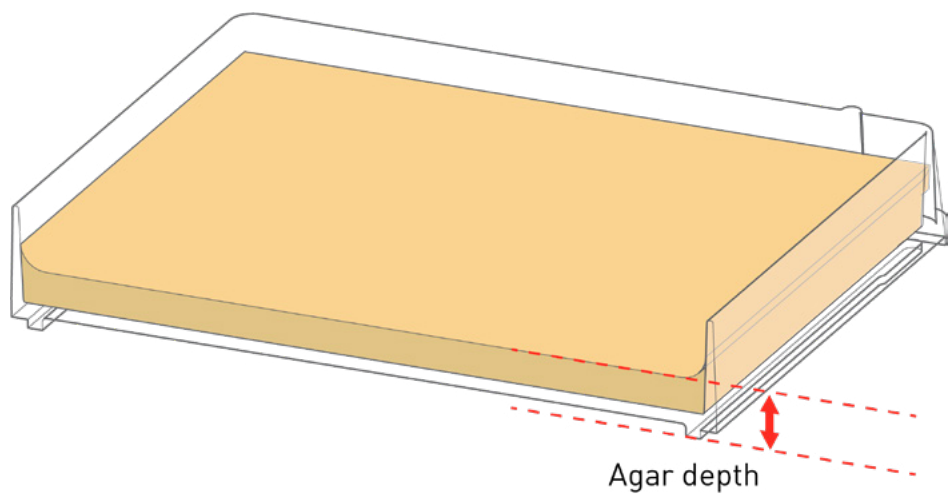



SINGER INSTRUMENTS

A RESPONSIBILITY TO SCIENCE!

Roadwater,
Watchet,
Somerset.
TA23 0RE
UK

+44 (0)1984 640226 (tel)
+44 (0)1984 641166 (fax)

contact@singerinstruments.com
singerinstruments.com



**SCAN TO VISIT WEBSITE FOR MORE
HELPFUL TIPS AND TUTORIALS!**



PHEN@BOOTH+